

Catalog Number: 100831, 100834, 191335, 195303

Lysozyme

CAS # : 9001-63-2

Synonym: Muramidase

Physical Appearance: White to off white lyophilized crystalline powder

Description: Lysozyme (muramidase) hydrolyzes preferentially the b-1,4 glucosidic linkages between N-acetylmuramic acid and N-acetylglucosamine which occur in the mucopeptide cell wall structure of certain microorganisms, such as *Micrococcus lysodeikticus*.

In humans, Osserman, et al.¹⁴ have indicated that lysozyme may be the mediator in the anti-tumor function of macrophages which, it has been shown, secrete the enzyme.⁸ There is evidence that cartilage lysozyme has a role in cartilage calcification.¹²

The enzyme has been used for lysing *E. coli* and *Streptomyces* for extraction purposes⁹ such as extracting group specific antigen.¹⁸ It would appear that lysozyme may act as a germinative agent of bacterial spores.^{2,7}

Molecular weight: 14,388 (Chicken¹¹); 17,000 (human).

Optimum pH: 9.2.⁵

Extinction coefficient: = 26.4.³

Isoelectric point: pH 11.0.¹

Inhibitors: The enzyme is inhibited by surface-active reagents such as dodecyl sulfate, alcohols and fatty acids.¹⁷ Imidazole and indole derivatives are inhibitors via formation of charge-transfer complexes.¹⁵

Stability: Solutions at pH 4-5 are stable for several weeks refrigerated and for days at ambient temperatures.

Assay: Mintz, et al.¹³ describe a sensitive fluorimetric assay.

Unit Definitions:

From Chicken: One unit will cause a decrease in A450 of 0.001 per minute at pH 6.24 and 25°C using *Micrococcus lysodeikticus* as substrate.

Shugar Unit:¹⁶ One unit is defined as the amount of enzyme that will digest powdered cells of *Micrococcus lysodeikticus*, causing a decrease in absorbancy of 0.001 per minute at 37°C, pH 7.0

Availability:

Catalog Number	Description	Size
100831	Lysozyme, from chicken egg white, 3X crystallized; Activity: 20,000 - 25,000 units/mg	100 mg 500 mg

	protein	1 g 5 g 10 g 25 g 100 g
100834	Lysozyme, from chicken egg white, 2X crystallized; Activity: > 9,000 units/mg protein	1 g 5 g 10 g
195303	Lysozyme, from chicken egg white, Type VI, 3X crystallized; Activity: ~60,000 units/mg protein	1 g 5 g 25 g
191335	Lysozyme, from human neutrophils; Activity: ~30,000 Shugar units/mg protein	100 ug

References:

1. Alderton, G., Ward, W., and Febold, H., "Isolation of Lysozyme from Egg White," *J. Biol. Chem.*, **v. 157**, 43 (1945).

2. Ando, Y., "The Effect of Lysozyme on Ionic Forms of Spores of *Clostridium perfringens* Type A," *J. Bacteriol.*, **v. 122**, 794 (1975).

3. Aune, K., and Tanford, C., "Thermodynamics of the Denaturation of Lysozyme by Guanidine Hydrochloride. I. Dependence on pH at 25°C.," *Biochem.*, **v. 8**, 4579 (1969).

4. Dahlquist, F., Borders, C., Jacobson, G., and Raftery, M., "The Stereospecificity of Human, Hen and Papaya Lysozymes," *Biochem.*, **v. 8**, 694 (1969).

5. Davies, A., Neuberger, A., and Wilson, B., "The Dependence of Lysozyme Activity on pH and Ionic Strength," *Biochim. Biophys. Acta*, **v. 178**, 294 (1969).

6. Dubois, T., Guillard, R., Prieels, J., and Perraudin, J., "Comparison between the Folding of Reduced Hen Egg White Lysozyme and that of Reduced Human Milk Lysozyme," *Biochem.*, **v. 21**, 6516 (1982).

7. Duncan, C., Labbe, R., and Reich, R., "Germination of Heat- and Alkali-Altered Spores of *Clostridium perfringens* Type A by Lysozyme and and Initiation Protein., " *J. Bacteriol.*, **v. 109**, 550 (1972).

8. Gordon, S., Todd, J., and Cohn, Z., "In Vitro Synthesis and Secretion of Lysozyme by Mononuclear Phagocytes," *J. Exp. Med.*, **v. 139**, 1228 (1974).

9. Haas, M., and Dowding, J., "Aminoglycoside-Modifying Enzymes," in *Methods in Enzymology*, **XLIII**, (Hash, J., ed.), Academic Press, NY, 621 (1965).

10. Hankiewicz, J., and Swierczek, E., "Lysozyme in Human Body Fluids," *Clin. Chim. Acta*, **v. 57**, 205 (1974).

11. Jollés, P., "Lysozyme," in *The Enzymes*, **IV**, 2nd Ed., (Boyer, P., Lardy, H., and Myrback, K., eds.), Academic Press, NY, 431 (1960).

12. Kuettner, K., Sorgente, N., Croxen, R., Howell, D., and Pita, J., "Lysozyme in Preosseous Cartilage. VII. Evidence for Physiological Role of Lysozyme in Normal Endochondral Calcificaion," *Biochim. Biophys. Acta*, **v. 372**, 335 (1974).

13. Mintz, G., Herbold, D., and Glaser, L., "A Fluorescent Assay for Bacterial Cell Wall Lytic Enzymes," *Anal. Biochem.*, **v. 66**, 272 (1975).

14. Osserman, E., Kockars, M., Halper, J., and Fischel, R., "Effects of Lysozyme on Normal and Transformed Cells," *Nature*, **v. 243**, 331 (1973).

15. Shinitzky, M., Katchalski, E., Grisaro, V., and Sharon, N., "Inhibition of Lysozyme by Imidazole and Indole Derivatives," *Arch. Biochem. Biophys.*, **v. 116**, 332 (1966).

16. Shugar, D., "Measurement of Lysozyme Activity and the Ultra Violet Inactivation of Lysozyme," *Biochim. Biophys. Acta*, **v. 8**, 302 (1952).

17. Smith, G., and Stoker, C., "Inhibition of Crystalline Lysozyme," *Arch. Biochem.*, **v. 21**, 383 (1949).

18. Watson, B., Moellering, R., and Kunz, L., "Identification of *Streptococci*: Use of Lysozyme and *Streptomyces albus* Filtrate in the Preparation of Extracts for Lancefield Grouping," *J. Clin. Microbiol.*, **v. 1**, 274 (1975).

