



Instruction Manual

E-Gel® 96 Editor Software

For Macintosh®

Version A
September 3, 2002
25-0613

Table of Contents

| | |
|---|-----------|
| Installation | 1 |
| Downloading and Installing E-Gel® 96 Editor for Macintosh® | 1 |
| Computer Requirements | 1 |
| QuickTime™ Requirement | 1 |
| Memory Settings..... | 2 |
| Introduction | 3 |
| Welcome to E-Gel® 96 Editor for Macintosh | 3 |
| Imaging Gels | 3 |
| Supported File Formats | 3 |
| Compatible Image Analysis Software | 3 |
| Cropping Images | 3 |
| Getting Started | 4 |
| Starting the Program | 4 |
| Workflow Overview | 4 |
| Opening and Saving Images | 5 |
| Opening Images..... | 5 |
| Saving Images | 5 |
| Adjusting the Image Display | 6 |
| Magnifying Images | 6 |
| Adjusting the Image Brightness | 6 |
| Aligning the Targets | 7 |
| Introduction | 7 |
| Rotating the Image | 7 |
| Aligning the Red Targets with the Gel Crosshairs | 8 |
| Completing the Alignment | 8 |
| Adjusting the Lane Boxes | 9 |
| Lane Boxes..... | 9 |
| Adjusting the Width or Horizontal Position of All Lane Boxes..... | 10 |
| Manually Repositioning One or More Lane Boxes..... | 11 |
| Undoing Lane Box Adjustments | 11 |
| Saving Alignment Settings..... | 11 |
| Completing the Adjustment | 11 |
| Printing, Copying, and Analyzing the Image | 12 |
| Displaying the Configured Lanes | 12 |
| Hiding and Displaying Lane Labels | 12 |
| Smoothing the Image | 13 |
| Copying the Entire Image | 13 |
| Copying Individual Lanes..... | 14 |
| Printing the Image..... | 14 |
| Further Analysis | 14 |
| Software License | 15 |
| Technical Service | 16 |
| World Wide Web | 16 |

Contact Us16
Limited Warranty17

Installation

Downloading and Installing E-Gel[®] 96 Editor for Macintosh[®]

The E-Gel[®] 96 Editor for Macintosh[®] installation file can be downloaded for free from the Invitrogen Web site.*

1. Go to the E-Gel[®] product page at www.invitrogen.com/egels and follow the links to the **E-Gel[®] 96 Editor Software download** page.
2. Follow the instructions on the page to download the Macintosh[®] installation file (**E-Gel 96 Editor.sit** or **E-Gel 96 Editor.hqx**) to your computer.
3. After downloading the file, double-click on it to unstuff the application. If your computer does not include the Stuffit[™] Expander utility, you can download a version for free from www.stuffit.com.
4. The **E-Gel 96 Editor** folder, containing the program and associated files, will be extracted into the same location as the .sit file. You can move this folder anywhere on your hard drive.

*If your computer is not connected to the Internet, you can download the file to another computer and copy it to a removable disk for installation. If you do not have Internet access, contact Technical Service (see page 16) to obtain a copy of the file.

Computer Requirements

The following are the minimum system requirements for running E-Gel[®] 96 Editor for Macintosh[®]:

Operating system: Mac OS 8.5 or higher (OS X w/ OS 9 emulation only)
Computer: G3 Power Mac or higher
Built-in memory: 128 MB or higher

QuickTime[™] Requirement


E-Gel[®] 96 Editor for Macintosh[®] requires that QuickTime[™] version 4.5 or later be installed on your Macintosh[®] to display images. QuickTime[™] can be downloaded for free from: <http://www.apple.com/quicktime/>

Installation, continued

Memory Settings

The default application memory setting for E-Gel® 96 Editor on the Macintosh® is designed for a typical image size of about 1 MB. If your images are larger than this, you may experience system memory problems using the application. You can increase your virtual memory allocation by following the steps below:

1. With the program closed, locate and select the **E-Gel 96 Editor** application icon on your Macintosh® hard drive (do not select a shortcut icon; select only the original program icon).
2. On the Macintosh® Desktop, go to the **File** menu and select **Get Info**, and then select **Memory**.
3. In the Memory settings dialog, set *both* the **Preferred Size** and **Minimum Size** to approximately 20 times the typical image size. For example, if your images are around 2 MB (2000 K), change both sizes to 40000 K.

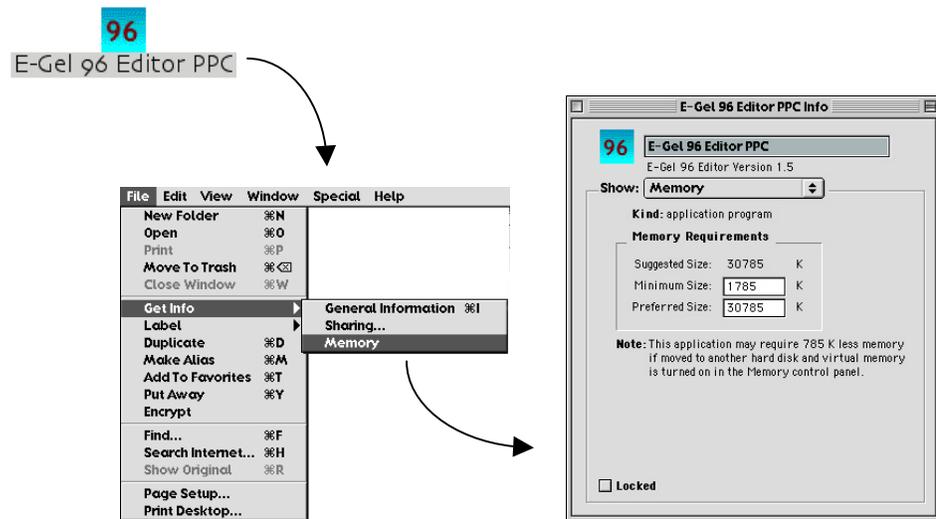

Note

- If the built-in memory on your Macintosh® is lower than 128 MB, increasing the virtual memory may not help with large images.
- A higher minimum virtual memory setting for E-Gel® 96 Editor may make it difficult for you to run other applications (which also require memory), particularly if your built-in system memory is 128 MB or less. Select the smallest memory setting necessary for the program to run smoothly.

Introduction

Welcome to E-Gel® 96 Editor for Macintosh

E-Gel® 96 Editor software for Macintosh® allows you to quickly reconfigure digital images of E-Gel® 96 gels for analysis and documentation. The staggered lanes in an E-Gel® 96 gel can be difficult to compare and analyze using standard 1-D gel analysis programs such as Bio-Rad's Quantity One, Phoretix 1D, or Kodak 1D software. E-Gel® 96 Editor reconfigures the staggered wells into a side-by-side format for easy comparison and analysis.

After scanning or capturing a digital image of an E-Gel® 96 gel, open the image in E-Gel® 96 Editor and follow the steps described in this manual to align and arrange the lanes in the image. Then save the reconfigured image for further analysis, print the image, and/or copy and paste selected lanes or the entire image into other applications for printing, saving, e-mailing, and/or publishing on the Web.

Imaging Gels

Use a flatbed scanner or digital camera to capture a digital image of your E-Gel® 96 gel. **Note:** When imaging, the gel should be properly aligned (i.e., not at an angle), and gel features should be clear and distinct in the image.



Important

The crosshairs ⊕ molded along the sides of the E-Gel® 96 cassette are essential for aligning the image in the software, and must be clear and distinct in the image. We recommend an image resolution of 150 dpi or higher.

Supported File Formats

After imaging, save the image in one of the file formats supported by E-Gel® 96 Editor for Macintosh® listed below:

- tiff
- bitmap
- gif
- jpeg
- png
- pict
- Quicktime

Note: If you will be analyzing the gel image using an analysis software program, be sure to save the image in a format supported by that program.

Compatible Image Analysis Software

Images edited by E-Gel® 96 Editor can be analyzed using many different biological image analysis programs, including Bio-Rad's Quantity One, Phoretix 1D, Kodak 1D, and UVP's LabWorks and GelWorks software. Consult your image analysis software documentation for the appropriate file formats.

Cropping Images

If there is uneven border space around the edges of the E-Gel® 96 image, E-Gel® 96 Editor may misalign the lanes. You may need to crop the border space using a standard image editing program such as Apple® Picture Viewer, QuickTime™ player, or Adobe® Photoshop®. Consult your image editing program documentation for detailed instructions.

Continued on next page

Getting Started

Starting the Program

To begin running E-Gel[®] 96 Editor, double-click on the application icon:



The graphical interface consists of a menu bar, an image window and toolbar, and a status bar.



Workflow Overview

E-Gel[®] 96 Editor software is designed to guide you through the following series of steps:

1. **Aligning the Image** (page 7). In the first step, you position red targets over the cross-hairs on the gel cassette to facilitate alignment.
2. **Adjusting Lanes** (page 9). In this step, you define the lanes on the gel for reconfiguration.
3. **Copying, Printing, and Analysis** (page 12). In the final step, the reconfigured image will be displayed on the screen, and you can proceed to copying, printing, and analysis.

In each step, use the **Next button** on the main toolbar to proceed to the next step, and the **Previous button** to return to the previous step:



Note

The image will not be reconfigured until you complete Step 2 above. When you complete Step 2, you should save the reconfigured image under a different name or in a different location if you want to preserve the original image.


Opening and Saving Images

Opening Images

To open an image in E-Gel[®] 96 Editor:

- Click on the **Open button** or select **Open** from the **File menu** (⌘+O) and select the image file using the **Open dialog** (see figure below); or
- Drag an image file directly into the open application window.

In the **Open dialog**, navigate to the folder containing your images and double-click on the image file name. The image will appear in the E-Gel[®] 96 Editor image window.

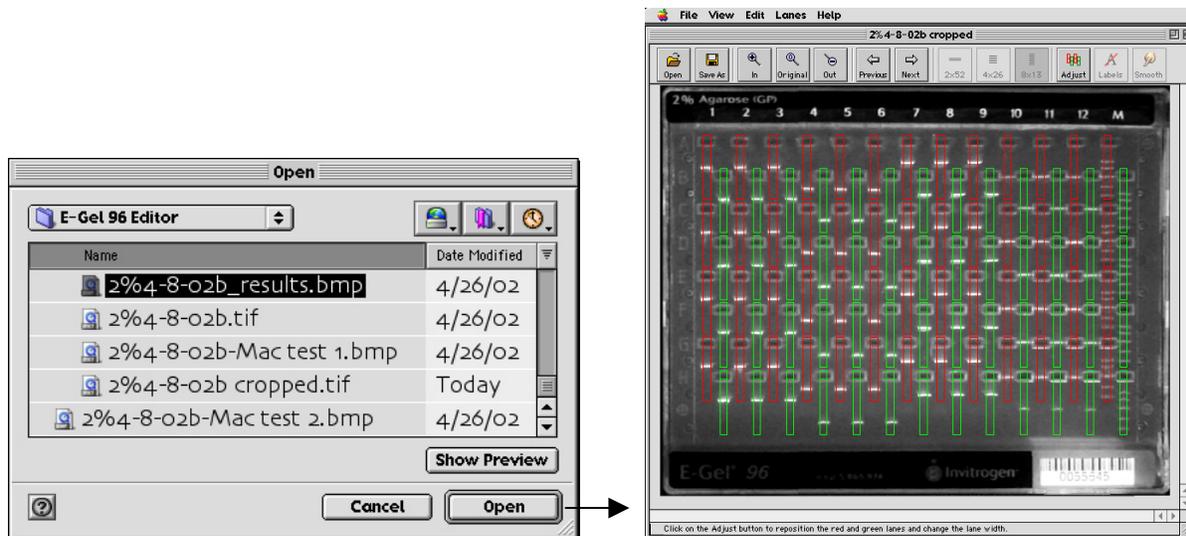

Note

- The E-Gel[®] 96 Editor window can contain only one image at a time; if you open another image, it will replace the current image. You will be prompted to save any changes to the current image.
- If you open an image that has already been reconfigured using E-Gel[®] 96 Editor, the **Previous** and **Next** buttons will be disabled, and only the options for reconfigured images will be available (changing the configuration, printing, copying, etc.). (See **Printing, Copying, and Analyzing the Image**, beginning on page 12.)

Saving Images

Click on the **Save As** button to open a dialog box in which you can save the image under the same name or a different name. This command is also on the **File menu**.

When saving images, note the following:

- E-Gel[®] 96 Editor will automatically append an underscore character (“_”) to the original image file name, to avoid overwriting the original file. To overwrite the original file, delete this character.
- E-Gel[®] 96 Editor automatically adds a file extension to the image file name, based on the file format (e.g., .jpg for jpeg images, .bmp for bitmaps).

To close an image without saving, select **Close** from the **File menu**.

Adjusting the Image Display

Magnifying Images

With the image open, click on the **Zoom In** button on the main toolbar to magnify the image, **Zoom Out** to shrink it, or **Original** to return the image to its actual size:




Adjusting the Image Brightness


If the image display is too bright or too dark, you can adjust it using the **Brightness** controls:

1. Go to the **View** menu and select **Brightness**.
2. In the **Brightness** dialog, drag the slider to make the image brighter or darker. To return to the original setting, click on the **Original** button.
3. When you are satisfied with the image brightness, click on **OK**. The image display will appear adjusted.


1. Image is too dark.

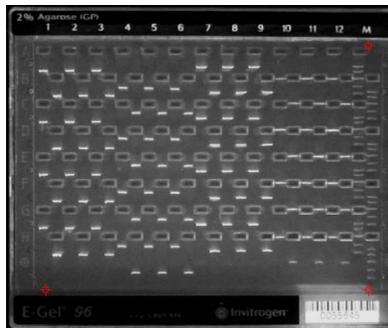

2. Open Brightness dialog, drag slider to lighten image, and click OK

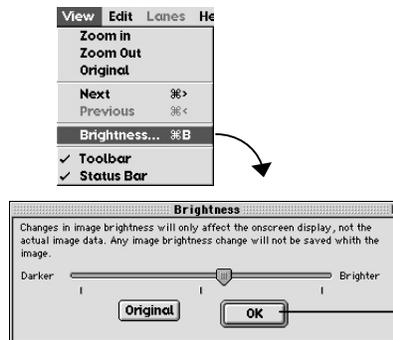

3. Image display is adjusted

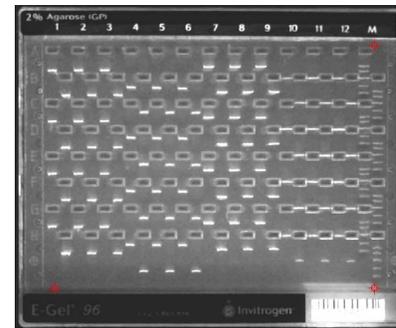

Aligning the Targets


Introduction

In this step, you prepare the image for alignment. First, make any necessary adjustments to the image orientation and brightness. Then position the three red targets displayed over the image directly over the three crosshairs ⊕ molded on the gel cassette.

Rotating the Image

If the E-Gel[®] 96 gel has been imaged at a 90° angle or upside-down, you can correct the orientation using the **Rotate** command.

Select **Rotate** from the **Edit menu**. The displayed image will rotate counterclockwise 90°. Select **Rotate** again to continue rotating the image.

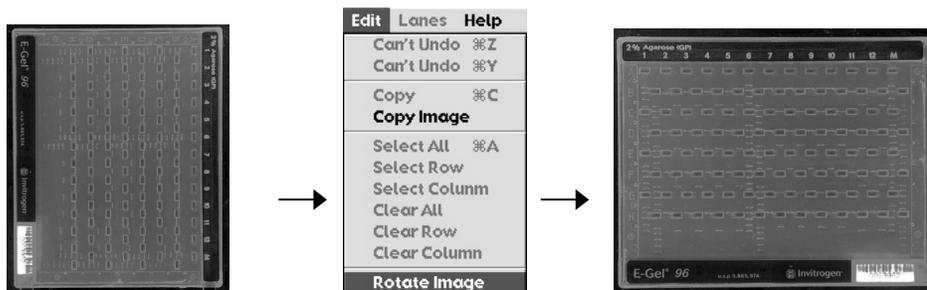

Note: This feature will change the image orientation; however, the image data (e.g., size and intensity of bands) will remain unchanged. The new orientation will be saved when you save the image.

Continued on next page

Aligning the Targets, continued


Aligning the Red Targets with the Gel Crosshairs

There are three red targets displayed over the image. To align the image, you must position these targets directly over the three crosshairs ⊕ molded on the lower left, upper right, and lower right corners of the gel cassette.


To do this, perform the following steps (see diagram below):

1. Locate a crosshair on one of the corners of the cassette. You can magnify the crosshair using the **Zoom In** tool to aid in target positioning.
2. Using your mouse:
 - **Drag** the nearest red target directly over the crosshair,
 - **Double-click** on the center of the crosshair to “jump” the target into position (you may need to fine-tune the position of the target after the jump), or
 - **Click on the target and use the arrow keys** to move the target over the crosshair.
3. Repeat this procedure for the remaining two crosshairs. Each red target should be directly aligned over a crosshair.


1. Magnify gel crosshair

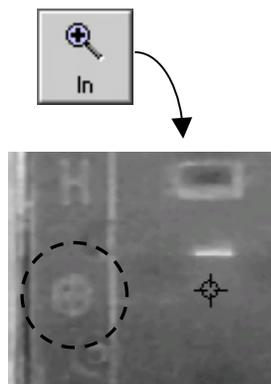

2. Drag, double-click, or use the arrow keys to move the target

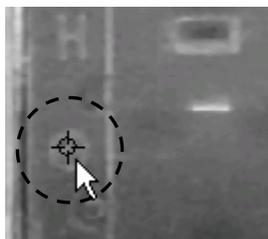

3. Repeat for all targets

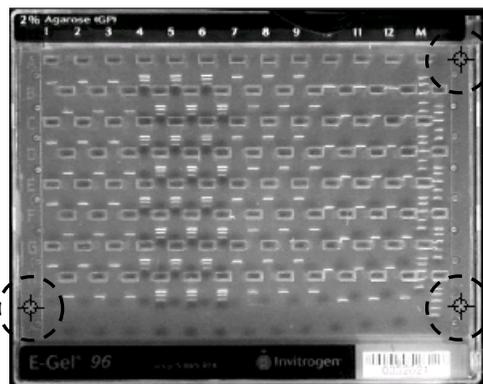

Important

The red targets must be positioned *directly* over the crosshairs for correct alignment.

Completing the Alignment

When the red targets are correctly positioned, click on the **Next button** to complete the alignment step.


Adjusting the Lane Boxes

Lane Boxes

When you complete the alignment step, red and green lane boxes will appear overlaid on the image (see the figure below).

These lane boxes indicate how the lanes will be cropped and positioned in the final configuration. In this step, you adjust the boxes so that they surround the individual lanes in the gel without overlapping adjacent lanes.

If the image is skewed or the lane boxes are grossly out of position, click on the **Previous button** to return to the alignment step.

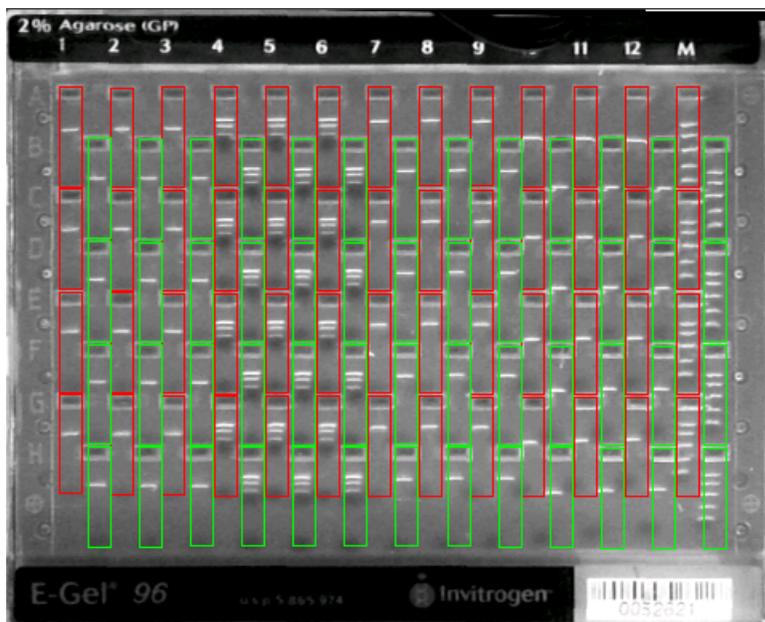

You can adjust all the lane boxes at the same time, or you can adjust one or more groups of boxes individually. See the following page for descriptions of both methods.



Important

If you analyze your gels using a 1-D gel image analysis software program, be careful to exclude the slopes of neighboring wells from the lane boxes. The software could misinterpret adjacent wells as bands.



Note

If you cannot successfully align the lane boxes, the image may contain border space around the edges, preventing proper alignment. Follow the steps below to remove this extra space:


1. Close the image in E-Gel[®] 96 Editor.
2. Open the image in an image-editing program such as Apple[®] Picture Viewer, QuickTime[™] player, or Adobe[®] Photoshop[®] and crop any extra space around the edges of the gel (see the program documentation for detailed instructions).
3. Re-open the image in E-Gel[®] 96 Editor and return to the alignment step.

Continued on next page

Adjusting the Lane Boxes, continued

Adjusting the Width or Horizontal Position of All Lane Boxes

To adjust the width and/or horizontal position of all of the lane boxes, click on the **Adjust** button or select the **Adjust** command from the **Lanes** menu to open the **Adjust Lanes** dialog box:

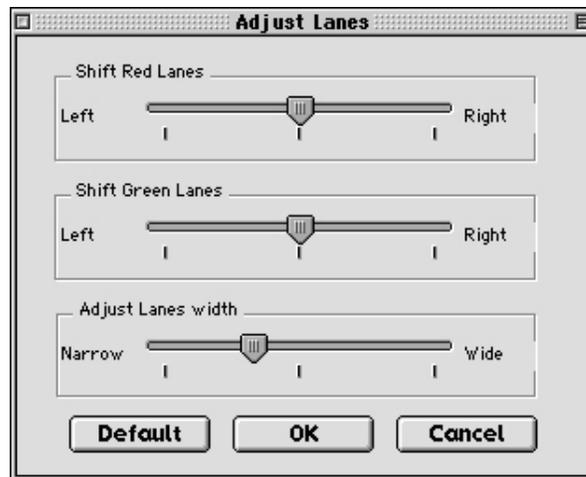

Use the **Shift Red Lanes** and **Shift Green Lanes** sliders to control the horizontal position of the lane overlays. Drag the sliders to move the red or green overlays left or right on the image.

Use the **Adjust Lane Width** slider to control the width of the lane overlays. Drag the slider to make the lanes wider or narrower.

The **Default** button resets the sliders to their default positions.

When the boxes are in the correct position, click on **OK** to close the dialog box.


Continued on next page

Adjusting the Lane Boxes, continued

Manually Repositioning One or More Lane Boxes

To adjust the horizontal and/or vertical position of one or more lane boxes:

1. Select the lane box or boxes:
 - To select lane boxes individually, click on them one at a time..
 - To select all the boxes in a column, select **Select Column** from the **Edit menu** and click on the column, or hold down the **Ctrl key**, click on the column, and select **Select Column** from the context menu.
 - To select all the boxes in a row, select **Select Row** from the **Edit menu** and click on the row, or hold down the **Ctrl key**, click on the row, and select **Select Row** from the context menu.
 - To select all the boxes in the image, select **Select All** from the **Edit menu** or hold down the **Ctrl key**, click on any lane in the image, and select **Select All** from the context menu.
2. Drag the lane box(es) to the new position, or use the arrow buttons on your keyboard to reposition the boxes.

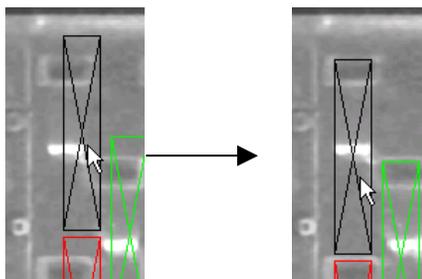

Undoing Lane Box Adjustments

To undo a manual lane box position change, select **Undo** from the **Edit menu** ($\text{⌘}+\text{Z}$). To redo the change, select **Redo** ($\text{⌘}+\text{Y}$).

Saving Alignment Settings

To save the aligned lane boxes for use on similar gel images:

1. Reconfigure the lane boxes as desired.
2. Select **Export Settings** from the **Lanes menu**.
3. In the **Export Settings dialog**, enter a name for the settings and save them in an appropriate location on your computer.

To recall saved settings on a new image:

1. With the image open and lane boxes displayed, select **Import Settings** from the **Lanes menu**.
2. In the dialog, select the settings file to open.
3. The lane boxes will be reconfigured based on the saved settings.


Completing the Adjustment

Confirm that the lane boxes completely surround the lanes in the gel without overlapping adjacent lanes. When you are satisfied with the position of the overlays, click on the **Next button** to complete the step.

Printing, Copying, and Analyzing the Image

Displaying the Configured Lanes

When you complete the adjust lanes step, the gel image will appear reconfigured in the application window:

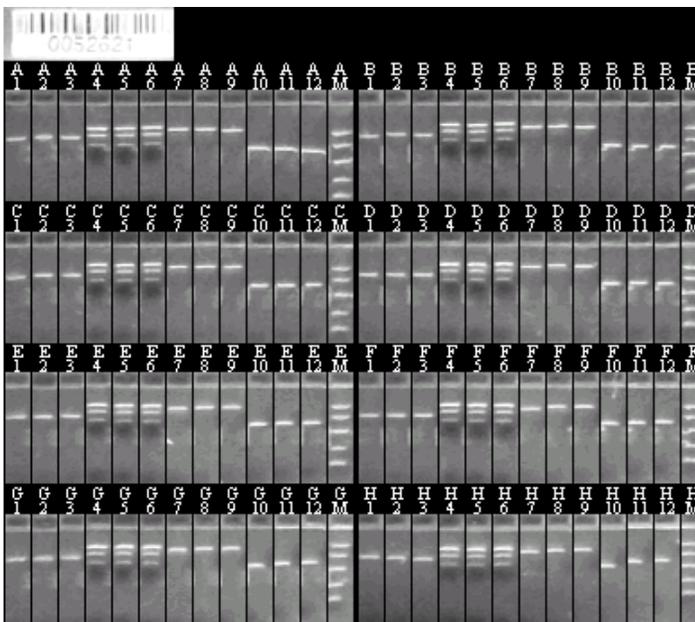


All of the lanes in each row will appear side-by-side for easy comparison. To change the configuration, click on the desired configuration button to rearrange the rows and columns in the image:



Hiding and Displaying Lane Labels

The labels marking each lane (A1, A2, A3, etc.) can be displayed or concealed for copying, printing, and publishing purposes. Click on the **Labels** button to hide the labels, and click on it again to display the labels.

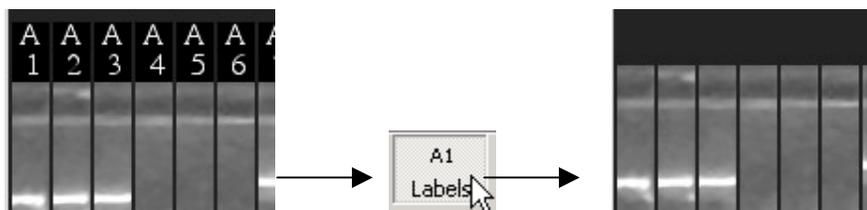

Continued on next page

Printing, Copying, and Analyzing the Image, continued

Smoothing the Image


Image smoothing may be desirable if the bands in the configured image appear jagged or speckled. Smoothing will remove minor imperfections in the image and create more uniform edges (see the example images below).

Note: Smoothing will permanently change the underlying image data; therefore, we strongly recommend that you save the smoothed image under a different file name for visual analysis, printing, or publishing only, and preserve the unsmoothed original for analysis using gel analysis software.

To perform smoothing:

1. Click on the **Smooth button**. You will be prompted to save the image under a different file name after smoothing is complete.
2. Click on **OK** to proceed.

Smoothing before and after:

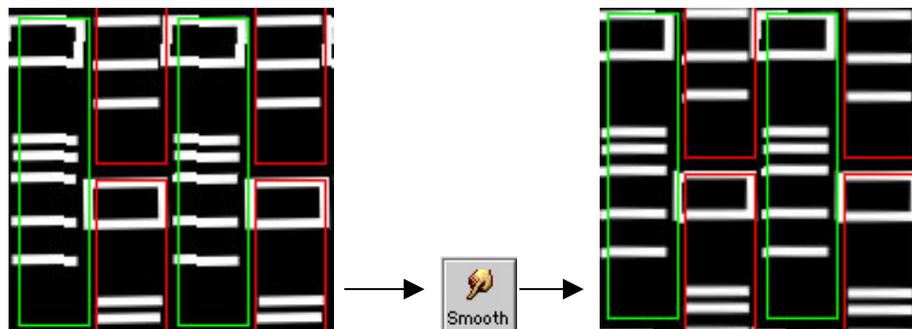

Note

Only the regions of the image defined by the lane boxes will be affected by the smoothing function. If you use the **Previous button** to go back to the unconfigured image, the areas defined by the lane boxes will appear smoothed, while the remaining areas will appear unchanged.

Copying the Entire Image

You can copy the entire image to the Macintosh® clipboard and then paste it into a document or other file. (To copy individual lanes, see below.)

To copy the entire image:

1. Select **Copy Image** from the **Edit menu** or hold down the **Ctrl key**, click on the image, and select the command from the context menu:



2. Open your word processing, spreadsheet, or other application and paste the image using the tools in that application.
-

Continued on next page

Printing, Copying, and Analyzing the Image, continued

Copying Individual Lanes


To select individual lanes in the image for copying, drag across the lanes or click on them one at a time. They will appear highlighted with a red border:



To deselect individual lanes, click on them again. To select all the lanes in the image, hold down the **Ctrl** key, click on the image, and select **Select All** from the context menu. To clear all selections, select **Clear All**.

When you are satisfied with your selections:

1. Select **Copy** from the **Edit** menu or context menu, or use the **⌘+C** key command. Only the selected lanes will be copied to the Macintosh® clipboard.
2. Paste the lanes into a word processing, spreadsheet, or other application for documentation and analysis. The lanes will be pasted as a block in alphanumerical order (regardless of the order of selection):

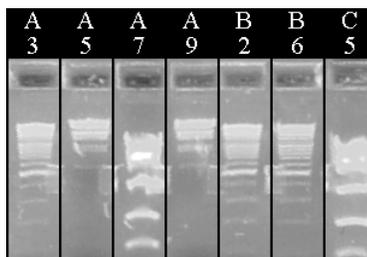

Printing the Image

To print the image, select **Print** from the **File** menu.

To change the printer and orientation settings, select **Page Setup** from the **File** menu.

Further Analysis

After you have reconfigured the image, click on the **Save As** button (**⌘+S**) to save it under a different name and open it in a 1-D gel analysis software program for further analysis.

Software License

Software License

By your use of E-Gel® 96 Editor Software, you accept and acknowledge the following terms and conditions:

- Invitrogen Corporation (“Invitrogen”) reserves all copyrights to E-Gel® 96 Editor Software.
- E-Gel® 96 Editor Software is free to customers who purchase E-Gel® 96 gels and E-Gel® 96 apparatus from Invitrogen.
- E-Gel® 96 Editor Software may be obtained at www.invitrogen.com/egels.
- E-Gel® 96 Editor Software will be used only with E-Gel® gels and E-Gel® 96 apparatus and will be used only for the internal research use of customers who purchase E-Gel® 96 gels and E-Gel® 96 apparatus from Invitrogen.
- E-Gel® 96 Editor Software will not be copied, transferred or redistributed.
- E-Gel® 96 Editor Software or its output will not be used for any COMMERCIAL use (including, but not limited to, selling, renting, leasing, or for use in the provision of a service or other commercial offering).

You agree not decompile, disassemble, reverse-engineer, or in any way modify the software or merge it into another program.

This software is provided “as is,” and any express or implied warranties including, but not limited to, the implied warranties of merchantability and fitness for a particular purpose are disclaimed. In no event shall invitrogen corporation be liable for any direct, indirect, incidental, special, exemplary, or consequential damages (including, but not limited to, procurement of substitute goods or services, loss of use, data, or profits, or business interruption) however caused and on any theory of liability, whether in contract, strict liability, or tort (including negligence or otherwise) arising in any way out of the use of this software, even if advised of the possibility of such damage. If any provision of this agreement is found to be legally unenforceable, then that provision will be deleted from this agreement, and the remainder of the agreement will continue in full force and effect.

Technical Service

World Wide Web



Visit the Invitrogen Web Resource using your World Wide Web browser. At the site, you can:

- Get the scoop on our hot new products and special product offers
- View and download vector maps and sequences
- Download manuals in Adobe® Acrobat® (PDF) format
- Explore our catalog with full color graphics
- Obtain citations for Invitrogen products
- Request catalog and product literature

Once connected to the Internet, launch your Web browser (Internet Explorer 5.0 or newer or Netscape 4.0 or newer), then enter the following location (or URL):

<http://www.invitrogen.com>

...and the program will connect directly. Click on underlined text or outlined graphics to explore. Don't forget to put a bookmark at our site for easy reference!

Contact Us

For more information or technical assistance, please call, write, fax, or email. Additional international offices are listed on our Web page (www.invitrogen.com).

Corporate Headquarters:

Invitrogen Corporation
1600 Faraday Avenue
Carlsbad, CA 92008 USA
Tel: 1 760 603 7200
Tel (Toll Free): 1 800 955 6288
Fax: 1 760 602 6500
E-mail:
tech_service@invitrogen.com

Japanese Headquarters:

Invitrogen Japan K.K.
Nihonbashi Hama-Cho Park Bldg. 4F
2-35-4, Hama-Cho, Nihonbashi
Tel: 81 3 3663 7972
Fax: 81 3 3663 8242
E-mail: jpinfo@invitrogen.com

European Headquarters:

Invitrogen Ltd
Inchinnan Business Park
3 Fountain Drive
Paisley PA4 9RF, UK
Tel: +44 (0) 141 814 6100
Tech Fax: +44 (0) 141 814 6117
E-mail:
eurotech@invitrogen.com

Continued on next page

Technical Service, Continued

Limited Warranty

Invitrogen is committed to providing our customers with high-quality goods and services. Our goal is to ensure that every customer is 100% satisfied with our products and our service. If you should have any questions or concerns about an Invitrogen product or service, please contact our Technical Service Representatives.

Invitrogen warrants that all of its products will perform according to the specifications stated on the certificate of analysis. The company will replace, free of charge, any product that does not meet those specifications. This warranty limits Invitrogen Corporation's liability only to the cost of the product. No warranty is granted for products beyond their listed expiration date. No warranty is applicable unless all product components are stored in accordance with instructions. Invitrogen reserves the right to select the method(s) used to analyze a product unless Invitrogen agrees to a specified method in writing prior to acceptance of the order.

Invitrogen makes every effort to ensure the accuracy of its publications, but realizes that the occasional typographical or other error is inevitable. Therefore Invitrogen makes no warranty of any kind regarding the contents of any publications or documentation. If you discover an error in any of our publications, please report it to our Technical Service Representatives.

Invitrogen assumes no responsibility or liability for any special, incidental, indirect or consequential loss or damage whatsoever. The above limited warranty is sole and exclusive. No other warranty is made, whether expressed or implied, including any warranty of merchantability or fitness for a particular purpose.

©2002 Invitrogen Corporation. All rights reserved.



United States Headquarters:

Invitrogen Corporation
1600 Faraday Avenue
Carlsbad, California 92008
Tel: 1 760 603 7200
Tel (Toll Free): 1 800 955 6288
Fax: 1 760 603 7229
Email: tech_service@invitrogen.com

European Headquarters:

Invitrogen Ltd
3 Fountain Drive
Inchinnan Business Park
Paisley PA4 9RF, UK
Tel (Free Phone Orders): 0800 269 210
Tel (General Enquiries): 0800 5345 5345
Fax: +44 (0) 141 814 6287
Email: eurotech@invitrogen.com

International Offices:

Argentina 5411 4556 0844
Australia 1 800 331 627
Austria 0800 20 1087
Belgium 0800 14894
Brazil 0800 11 0575
Canada 800 263 6236
China 10 6849 2578
Denmark 80 30 17 40

France 0800 23 20 79
Germany 0800 083 0902
Hong Kong 2407 8450
India 11 577 3282
Italy 02 98 22 201
Japan 03 3663 7974
The Netherlands 0800 099 3310
New Zealand 0800 600 200
Norway 00800 5456 5456

Spain & Portugal 900 181 461
Sweden 020 26 34 52
Switzerland 0800 848 800
Taiwan 2 2651 6156
UK 0800 838 380

For other countries see our website

www.invitrogen.com

