Cell Explorer[™] Live Cell Tracking Kit *Green Fluorescence*

Ordering Information	Storage Conditions	Instrument Platform
Product Number: 22621 (2 plates)	Keep in freezer Protect from moisture and light	Fluorescence microscope Flow cytometer

Introduction

Our Cell ExplorerTM Live Cell Tracking Kits are a set of tools that provide well-retained cell-tracing reagents to label and track cells for investigations of cellular functions in a wide variety of colors. The effective labeling of cells provides a powerful method for studying cellular events in a spatial and temporal context. This particular kit is designed to label live cells in green fluorescence for the studies that require the fluorescent tag molecules retained inside cells for a relatively longer time. The Track ItTM Green used in the kit is a non-fluorescent dye that carries a cell-retaining moiety. The dye becomes strongly fluorescent upon entering into live cells, and trapped inside cells to give a stable fluorescence signal. The adduct that formed in labeled cells is retained by the cells throughout development and meiosis, and is inherited by daughter cells after cell division.

This kit can be readily adapted for many different types of fluorescence platforms such as microplate assays, and flow cytometry analysis. For the long term tracking of cell proliferation or cell division studies, Track ItTM Green has the similar features as traditional cell proliferation reagent CFSE, but eliminates the dye efflux drawback associated with CFSE. The CFSE-stained cells show continued efflux of the dye molecules after 24 hours and the fluorescence intensity of daughter generations reduces by >10 times from first generation. Another advantage of Track ItTM Green is that it is compatible with cell culture medium in the staining cells prior to imaging or flow cytometric analysis, and those features make this kit more convenient and less hands-on operations than CFSE staining. The green fluorescent indicator used in the kit has Ex/Em = 490/520 nm, compatible with the FITC filter set that is installed with almost every major fluorescence instrument and flow cytometer.

Key Features		
Sensitive:	Detect at least 10 passages of Jurkat cells, 3 passages more than CFSE.	
Robust:	Uniform fluorescence distribution from generation to generation.	
<i>Rapid:</i> No efflux effect at first passage as CFSE does.		
Convenient:	No need to remove cell culture medium in the staining cells.	

Kit Components

Components	Amount
Component A: Track It [™] Green	1 vial
Component B: Assay Buffer	1 bottle (25 mL)
Component C: DMSO	1 vial (100 μL)

Assay Protocol

Brief Summary

Prepare samples → Add Track ItTM Green working solution → Stain the cells at RT for 15 to 30 minutes → Wash the cells → Examine the specimen at Ex/Em = 490/520 nm

Note: Thaw all the components at room temperature before opening.

1. Prepare Cells:

- 1.1 <u>For adherent cells</u>: Plate cells overnight in growth medium at 10,000 to 40,000 cells/well/90 μL for 96-well plates or 2,500 to 10,000 cells/well/20 μL for 384-well plates.
- 1.2 For non-adherent cells: Centrifuge the cells from the culture medium and then suspend the cell pellets in culture medium at 50,000-100,000 cells/well/90 μL for 96-well poly-D lysine plates or 10,000-25,000 cells/well/20 μL for 384-well poly-D lysine plates. Centrifuge the plates at 800 rpm for 2 minutes with brake off prior to the experiments.

Note 1: For flow cytometry experiment, prepare cells in 0.5 mL warm medium or buffer of your choice at a density of 5×10^5 to 1×10^6 cells/mL.

Note 2: Each cell line should be evaluated on an individual basis to determine the optimal cell density.

2. Prepare Track ItTM Green stain solution:

- 2.1 <u>Prepare 1000X Track ItTM Green stock solution</u>: Add100 µL of DMSO (Component C) into the vial of Track ItTM Green (Component A) and mix well. Note: The unused portion of 1000X Track ItTM stock solution should be stored at -20°C. Avoid repeated freeze/thaw cycles.
- 2.2 Prepare Track It[™] Green working solution: Dilute 1000X Track It[™] Green stock solution (from Step 2.1) into Assay Buffer (Component B) at 1:1000 ratio to make Track It[™] Green_working solution. Note: The final concentration of the Track It[™] Green should be empirically determined for different cell types and/or experimental conditions. In general, long-term staining (more than about 3 days) or the use of rapidly dividing cells will require 1: 500 dilution to double the dye concentration. Dye at a lower concentration up to 1:2000 dilution may be needed for shorter experiments, such as viability assays. To maintain normal cellular physiology and reduce potential artifacts, the concentration of the dye should be kept as low as possible.

3. Stain the cells:

- 3.1 Add equal volume of Track It[™] Green working solution (from Step 2.2) into the cell wells. For example, for 96-well plate, add 100 µL/well of Track It[™] Green working solution into the cells.
- 3.2 Incubate the cells in a 37 $^{\circ}$ C, 5% CO₂ incubator for 15 to 30 minutes.
- 3.3 Wash cells with HHBS or an appropriate buffer for 3 times. Note: Alternatively, fix the cells at this point. Store the fixed cells at 4 °C, and image the cells later.
- 3.4 Image the cells using a fluorescence microscope with FITC filters (Ex/Em = 490/520 nm). Or monitor the fluorescence intensity with a flow cytometer using the FL1 channel (Ex/Em = 490/525 nm), gate on the cells of interest, excluding debris.



Figure 1. Image of U2OS cells stained with Cell Explorer[™] Live Cell Tracking Kit in a Costar black 96-well plate.



Figure 2. Cell proliferation assay with Cell Explorer[™] Live Cell Tracking Kits. Jurkat cells were stained with Track It[™] Green or CFSE (0.5uM) on Day 0. The sample was allowed to proliferate for 9 days. Fluorescence intensity was measured with FACSCalibur (BD, San Jose, CA) flow cytometer in FL1 channel.

References

- 1. Wolff M, Wiedenmann J, Nienhaus GU, Valler M, Heilker R. (2006) Novel fluorescent proteins for highcontent screening. Drug Discov Today, 11, 1054.
- 2. Lee S, Howell BJ. (2006) High-content screening: emerging hardware and software technologies. Methods Enzymol, 414, 468.
- 3. Haasen D, Schnapp A, Valler MJ, Heilker R. (2006) G protein-coupled receptor internalization assays in the high-content screening format. Methods Enzymol, 414, 121.
- 4. Hudson CC, Oakley RH, Sjaastad MD, Loomis CR. (2006) High-content screening of known G proteincoupled receptors by arrestin translocation. Methods Enzymol, 414, 63.
- 5. Martinez ED, Dull AB, Beutler JA, Hager GL. (2006) High-content fluorescence-based screening for epigenetic modulators. Methods Enzymol, 414, 21.
- 6. Giuliano KA. (2007) Optimizing the integration of immunoreagents and fluorescent probes for multiplexed high content screening assays. Methods Mol Biol, 356, 189.
- 7. Gough AH, Johnston PA. (2007) Requirements, features, and performance of high content screening platforms. Methods Mol Biol, 356, 41.
- 8. Hoffman AF, Garippa RJ. (2007) A pharmaceutical company user's perspective on the potential of high content screening in drug discovery. Methods Mol Biol, 356, 19.
- 9. Taylor DL. (2007) Past, present, and future of high content screening and the field of cellomics. Methods Mol Biol, 356, 3.
- Wolff M, Haasen D, Merk S, Kroner M, Maier U, Bordel S, Wiedenmann J, Nienhaus GU, Valler M, Heilker R. (2006) Automated high content screening for phosphoinositide 3 kinase inhibition using an AKT 1 redistribution assay. Comb Chem High Throughput Screen, 9, 339.

Warning: This kit is only sold to end users. Neither resale nor transfer to a third party is allowed without written permission from AAT Bioquest®. Chemical analysis of the kit components is strictly prohibited. Please call us at 408-733-1055 or e-mail us at info@aatbio.com if you have any questions.