

ELISA diluent

- Product code:** 3652-D2
- Contents:** Protein-containing buffer, ready-to-use, with 0.15% Kathon CG as preservative.
- Applications:** To prevent interference by heterophilic antibodies in serum and plasma samples in ELISA. The diluent is recommended for use with human GM-CSF, Granzyme B, IFN- α , IFN- γ , IgE, IL-1 β , IL-2, IL-4, IL-5, IL-6, IL-10, IL-12, IL-13, IL-17A, IL-21, IL-23, IL-31, Perforin, Latent TGF- β 1, TNF- α , monkey IFN- γ , IL-2 and IL-17A ELISA tests. Please note that the ELISA diluent is optimized for use with antibodies from Mabtech.
- Instructions:** The ELISA diluent should be used for dilution of samples, cytokine standard and detection antibody. All samples should be diluted at least 2x in ELISA diluent. Do not use undiluted serum/plasma samples in the assay. Sample dilution can be made either in separate tubes or directly in the plate. If sample dilution is made directly in the plate, ELISA diluent should be added prior to addition of the sample.
- Important!** Do not use the ELISA diluent for dilution of Streptavidin conjugates. We recommend the use of an incubation buffer e.g. PBS with 0.05% Tween 20 containing 0.1% BSA for dilution of conjugates.
- Storage:** Shipped at ambient temperature. On arrival store product at 4-8°C. Adjust to room temperature before use.



Note; for research use only.

Mabtech shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.



MABTECH AB
Box 1233
SE-131 28 Nacka Strand
Sweden
Tel +46 8 716 27 00
Fax +46 8 716 27 01
mabtech@mabtech.com
www.mabtech.com

Mabtech, Inc.
M.E.B. Suite 220
3814 West Street
Mariemont, OH 45227
USA
Toll Free: +1 866 ELI-SPOT
Tel: +1 513 871 4500
Fax: +1 513 871 7353
mabtech.usa@mabtech.com

MABTECH AUSTRALIA
Pty Ltd
resolvingIMAGES
Unit 22
Thomastown Victoria 3074
Australia
Tel: +61 3 9466 4007
Fax: +61 3 9466 4003
mabtech.au@mabtech.com

Mabtech AB
Büro Deutschland
GERMANY
Tel: +49 40 4135 7935
Fax: +49 40 4135 7945
mabtech.de@mabtech.com

Mabtech AB
Bureau de liaison France
BP 255, 1300 route des
Crêtes
06905 Sophia Antipolis
FRANCE
Tel: +33 4 92 38 80 70
Fax: +33 4 92 38 80 71
mabtech.fr@mabtech.com