

## Purified anti-human IL-23 (p19)

**Catalog #/** 511201 / 50 µg

**Size:** 511202 / 500 µg

**Clone:** HLT2736

**Isotype:** Mouse IgG1, κ

**Immunogen:** Human recombinant p40 and p19 with elastin linker

**Reactivity:** Human, IL-23 p19 monomer, does not cross-react with mouse IL-23

**Preparation:** The antibody was purified by affinity chromatography.

**Formulation:** Phosphate-buffered solution, pH 7.2, containing 0.09% sodium azide.


**Storage:** The antibody solution should be stored undiluted at 4 °C.

## Applications


**Applications:** WB, IHC(Paraffin) (FPE using catalyzed signal amplification), IP

**Application Notes:** The purified HLT2736 antibody has been shown to be useful for Western blotting, immunohistochemistry using formalin-fixed, paraffin embedded tissues with catalyzed signal amplification), and immunoprecipitation.


**Recommended Usage:** Each lot of this antibody is quality control tested by ELISA assay. For ELISA capture applications, a concentration range of 2-8 µg/ml is recommended. To obtain a linear standard curve, serial dilutions of IL-23 recombinant protein ranging from 4000 to 30 pg/ml are recommended for each ELISA plate. It is recommended that the reagent be titrated for optimal performance for each application.

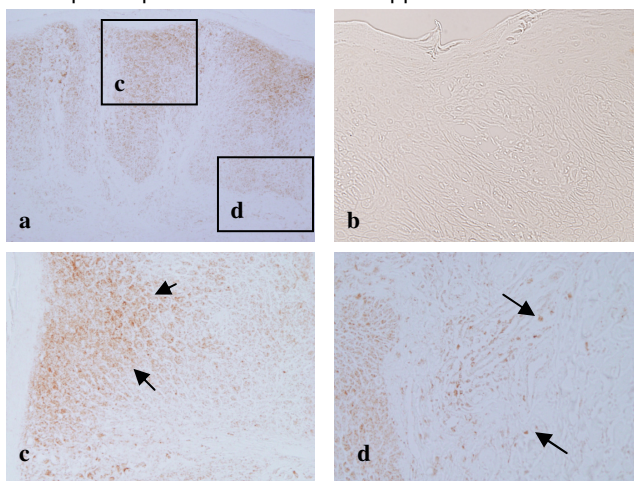

**Figure 1 Immunohistochemical localization of IL-23 expression in psoriatic lesions.** Paraffin embedded psoriatic lesion sections were stained with purified anti-human IL-23 (clone HLT2736) (**a, c & d**) and purified mouse IgG1,κ (clone MOPC-21) (**b**) as indicated. The brown color represents diaminobenzidine (DAB) deposition at the site of antibody binding. Low (×100, **a**) and high (×200, **b, c & d**) magnification are presented.

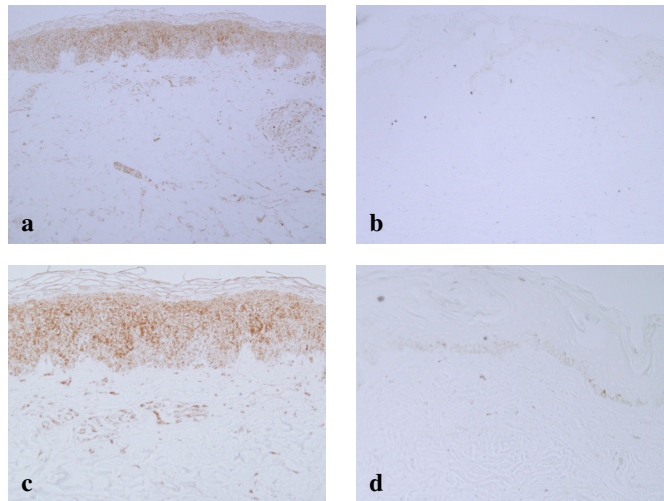

**Figure 2 IL-23 highly expressed in the uninvolved skin from psoriasis patients, but not in the control normal skin.** Paraffin embedded sections were stained with purified anti-human IL-23 (clone HLT2736) on the uninvolved skin from one psoriatic patient (**a & c**) and the control normal skin (**b & d**). Low ( $\times 100$ , **a & b**) and high ( $\times 200$ , **c & d**) magnification are presented.

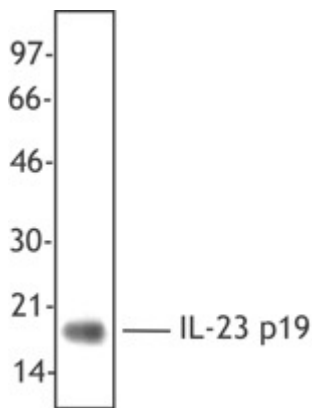

**Figure 3:** Linker-free, recombinant human IL-23 was resolved by electrophoresis, transferred to nitrocellulose and probed with the anti-IL23 p19 specific antibody HLT2736 antibody. Proteins were visualized using goat anti-mouse IgG conjugated to HRP and a chemiluminescence detection system. The HLT2736 recognizes the p19-specific subunit of human IL-23. This antibody will also recognize human IL-23 (p19 +p40 subunits) in recombinant IL-23 containing a linker moiety.

## Antigen Information

**Other Names:** p19p40, JKA3 induced upon T cell activation

**Structure:** Member of the IL-12 cytokine family, heterodimeric cytokine composed of p40 subunit shared with IL-12 and a specific p19 subunit

**Distribution:** Secreted by activated dendritic cells and macrophages

**Function:** Enhances IFN- $\gamma$  secretion by memory (CD45RO) T cells in an IL-2 dependent manner; may induce unique Th subset (designated Th<sub>IL-17</sub>) secreting IL-17, IL-6, TNF and low levels of IFN- $\gamma$  in EAE model; shown to promote immunity to mycobacteria

**Ligands/Receptors:** IL-23 receptor ( $\beta 1$  subunit of IL-12 receptor and IL-23-specific subunit)

**Regulation:** Upregulated in viral infections such as HSV, increased in psoriatic lesions in monocytes and mature dendritic cells

**Description:** The HLT2736 clone recognizes human IL-23, a heterodimeric cytokine composed of the p40 subunit shared with IL-12 and a specific p19 subunit. This antibody recognizes heterodimeric IL-23, but does not react with the p40 subunit alone or recombinant IL-12 when used as a capture antibody in ELISA. IL-23 is secreted by activated dendritic cells and macrophages and has been shown to enhance IFN- $\gamma$  secretion by memory (CD45RO) T cells in an IL-2 dependent manner. IL-23 may also induce unique Th subset (designated Th<sub>IL-17</sub>) that secretes the cytokines IL-17, IL-6, TNF and low levels of IFN- $\gamma$ . IL-23 has been shown to promote immunity to mycobacteria and has been shown to be upregulated in response to some viral infections and in psoriatic lesions. IL-23 binds to the IL-23 receptor comprised of a  $\beta$ 1 subunit shared with the IL-12 receptor and IL-23-specific subunit. This antibody has been shown to be useful for Western blotting, immunohistochemistry using formalin-fixed, paraffin embedded tissues with catalyzed signal amplification), and ELISA as a capture antibody.

**Antigen References:**

1. Parham, C., *et al.*, 2002. *J. Immunol.* 168:5699.
2. Verreck, F. A., *et al.*, 2004. *Proc. Natl. Acad. Sci. USA* 101:4560.
3. Lee, E., *et al.*, 2004. *J. Exp. Med.* 199:125.
4. Oppmann, B. *et al.*, 2000. *Immunity* 13:715.

---

## Protocol for IL-23 staining on paraffin embedded sections

**Reagents:**

1. Catalyzed Signal Amplification (CSA) System from DakoCytomation for use with mouse primary antibodies (Code K1500)
2. The CSA Ancillary System (Code K1499, DakoCytomation):
  - Target Retrieval Solution
  - Biotin Blocking Systems
  - TRIS-Buffered Saline with Tween 20 (TBST)(S-3306)
3. Purified anti-human IL-23 (clone HLT2736)
4. Purified mouse IgG1,k (clone MOPC-21)

**Staining Procedure:**

1. De-paraffinization of tissue sections: Prior to staining, deparaffinize tissue slides with xylene and gradient alcohols to remove embedding media and then re-hydrate.
2. Immerse antigen retrieval: tissue sections with Target Retrieval Solution (included in CSA Ancillary System) and microwave 1 min with high power and 9 min with medium power, then cool at room temperature for 20 min.
3. Blocking of endogenous avidin-binding activity (endogenous biotin): Sequentially apply avidin and biotin to tissue sections with Biotin blocking system (included in CSA Ancillary System).

4. Staining protocol:

- 1) Peroxidase block: Tap off excess liquid and wipe around specimen to remove any remaining liquid and to keep reagent within the prescribed area.

Apply enough hydrogen peroxide from bottle 1 (CSA System) to cover specimen, and incubate 5 ( $\pm$ 1) min.

Rinse gently with TBST buffer solution, and place in fresh bath for 3 min.

- 2) Protein block: Tap off excess liquid and wipe slide as before. Apply enough drops from bottle 2 (protein block, CSA System) to cover specimen. Incubate 5 ( $\pm$ 1) min. Do not rinse off protein block.

- 3) Primary antibody or negative control reagent: Tap off excess liquid and wipe slide as before.

Apply enough IL-23 antibody (20 µg/ml) or negative control IgG,κ to cover specimen. Incubate 15 (±1) min.  
Rinse gently with TBST buffer and place in up to three fresh TBST buffer baths for 3-5 minutes each.

4) Link antibody:

Tap off excess liquid and wipe slide as before.

Apply enough drops from bottle 3 (link antibody, CSA System) to cover specimens.

Incubate 15 (±1) min.

Rinse gently with TBST buffer and place in up to three fresh TBST buffer baths for 3-5 minutes each.

5) Streptavidin-biotin complex

Tap off excess liquid and wipe slide as before.

Apply enough drops of the prepared Streptavidin-Biotin Complex (CSA System) to cover specimens.

Incubate 15 (±1) min.

Rinse gently with TBST buffer and place in up to three fresh TBST buffer baths for 3-5 minutes each.

6) Amplification reagent

Tap off excess liquid and wipe slide as before.

Apply enough drops from Bottle 8 (Amplification reagent, CSA System) to cover specimens.

Incubate 15 (±1) min.

Rinse gently with TBST buffer and place in up to three fresh TBST buffer baths for 3-5 minutes each.

7) Streptavidin-peroxidase

Tap off excess liquid and wipe slide as before.

Apply enough drops from Bottle 9 (Streptavidin-Peroxidase, CSA System) to cover specimens.

Incubate 15 (±1) min.

Rinse gently with TBST buffer and place in up to three fresh TBS buffer baths for 3-5 minutes each.

8) Substrate-Chromogen solution

Tap off excess liquid and wipe slide as before.

Apply enough of the prepared substrate-Chromogen solution (CSA System) to cover specimens.

Incubate 15 (±1) min.

Rinse gently with TBST buffer and place in up to three fresh TBST buffer baths for 3-5 minutes each.

9) Mounting

Dehydrate specimens with gradient alcohol and xylene and mount with permanent mounting medium.



For research use only. Not for diagnostic use. Not for resale. Biolegend will not be held responsible for patent infringement or other violations that may occur with the use of our products.

BioLegend 11080 Roselle St., San Diego, CA 92121 [www.biolegend.com](http://www.biolegend.com)  
Toll-Free Phone: 1-877-Bio-Legend (246-5343) Phone: (858) 455-9588 Fax: (858) 455-9587