

# Magnetic Stand-96

Catalog Number AM10027

Pub. No. 4386455 Rev. C

Description	Quantity
Magnetic Stand-96	1

 **WARNING!** For every chemical, read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## Product description

The Magnetic Stand-96 is designed for paramagnetic bead precipitation from standard 96-well U-bottom microplates and 0.2-mL PCR plates with no additional accessories. The stand has a clear polyacryl base in standard SBS 96-well microplate footprint. The use of 24 extremely strong permanent magnets guarantees easy and fast magnetic separation in as little as 30 seconds. It sits flat on a lab bench as well as in a microplate adapter on robotic workstations.

## Using the Magnetic Stand-96

The Magnetic Stand-96 is ideal for use with MagMAX™-96 magnetic bead-based RNA isolation kits, MagMAX™ for Stabilized Blood Tubes RNA Isolation Kits, MagMAX™ FFPE Nucleic Acid Isolation Kits, and the MessageAmp™ Premier RNA Amplification Kit. For more information on these kits, visit their web catalog pages at [www.lifetechnologies.com](http://www.lifetechnologies.com).

**IMPORTANT!** We recommend cleaning the Magnetic Stand-96 with deionized water only; alcohol or alcohol-based cleaning agents may cause cracking of the stand material.

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions). If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

**For Research Use Only. Not for human or animal therapeutic or diagnostic use.**

The information in this guide is subject to change without notice.

**DISCLAIMER**

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

**NOTICE TO PURCHASER: LIMITED USE LABEL LICENSE: Research Use Only**

The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact [outlicensing@lifetech.com](mailto:outlicensing@lifetech.com) or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

**TRADEMARKS**

The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

© 2012 Life Technologies Corporation. All rights reserved.

**For support visit** [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support) or email [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

[www.lifetechnologies.com](http://www.lifetechnologies.com)