



Princeton Business Park, 5 Crescent Ave., P O Box 275
Rocky Hill, NJ 08553 USA
Tele: 609.497.0253 Fax: 609.497.0321
www.peprotech.com Email: info@peprotech.com

Murine IFN- γ
ELISA Development Kit
900-K98 Lot# 0212098
Expiration one year from date of receipt

Description: Murine IFN- γ ELISA development kit contains the key components required for the quantitative measurement of natural and/or recombinant mIFN- γ in a sandwich ELISA format within the range of 15–2000pg/ml. Using the ELISA protocol described below, the recommended microplates, reagents and solutions, the components supplied in this kit are sufficient to assay mIFN- γ in approximately 1000 ELISA plate wells.

RECONSTITUTION & STORAGE

Capture Antibody: 100 μ g of antigen-affinity purified rabbit anti-mIFN γ + 2.5mg D-mannitol. Centrifuge vial prior to opening. Reconstitute in 1.0ml sterile water for a concentration of 100 μ g/ml.

Detection Antibody: 25 μ g of biotinylated antigen-affinity purified rabbit anti-mIFN- γ + 2.5mg D-mannitol. Centrifuge vial prior to opening. Reconstitute in 0.25ml sterile water for a concentration of 100 μ g/ml.

Murine IFN- γ : 1 μ g of recombinant mIFN- γ + 2.2mg BSA + 11.0mg D-mannitol. Centrifuge vial prior to opening. Reconstitute in 1ml sterile water for a concentration of 1 μ g/ml.

Note: The reconstituted components are stable for 2 weeks when stored at 2–8°C. Components that have been reconstituted and aliquoted can be stored at –20°C for up to 6 months.

Avidin-HRP Conjugate: 60 μ l vial. Upon receipt, avidin-HRP conjugate should be aliquoted into ten 6 μ l vials and stored at \leq –20°C. Aliquots stored frozen at \leq –20°C are stable for up to 2 years from date of receipt. Avoid more than one freeze-thaw cycle. **Avidin should be used in conjunction with ABTS only.**

**RECOMMENDED MATERIALS (or purchase
PeproTech's ELISA Buffer Kit: Cat. # 900-K00)**

ELISA microplates (Nunc MaxiSorp Prod. # 439454, or Corning Prod # 3590);
Tween-20 (Sigma Cat. # P-7949);
BSA (Sigma Cat # A-7030);
ABTS Liquid Substrate Solution (Sigma Cat. # A3219);
Dulbecco's PBS [10x] (Gibco BRL Cat. # 14200-075).

RECOMMENDED SOLUTIONS

All solutions should be at ambient temperature prior to use.

PBS: dilute 10xPBS to 1xPBS, pH 7.20 in sterile water.

Wash Buffer: 0.05% Tween-20 in PBS

Block Buffer: 1% BSA in PBS *

Diluent: 0.05% Tween-20, 0.1% BSA in PBS *

*** Sterile filter and store at 4°C for up to 1 week.**

PLATE PREPARATION

1. Dilute capture antibody with PBS to a concentration of 1.0 μ g/ml. Immediately, add 100 μ l to each ELISA plate well. Seal the plate and incubate overnight at room temperature.
2. Aspirate the wells to remove liquid and wash the plate 4 times using 300 μ l of wash buffer per well.

After the last wash invert plate to remove residual buffer and blot on paper towel.

3. Add 300 μ l block buffer to each well. Incubate for at least 1 hour at room temperature.
4. Aspirate and wash plate 4 times.

ELISA PROTOCOL

Standard/Sample: Dilute standard from 2ng/ml to zero in diluent. Immediately add 100 μ l of standard or sample to each well in triplicate. Incubate at room temperature for at least 2 hours.

Detection: Aspirate and wash plate 4 times. Dilute detection antibody in diluent to a concentration of 0.25 μ g/ml. Add 100 μ l per well. Incubate at room temperature for 2 hours.

Avidin Peroxidase: Aspirate and wash plate 4 times. Dilute one 5.5 μ l aliquot of Avidin Peroxidase 1:2000 in diluent for total volume of 11ml. Add 100 μ l per well. Incubate 30 minutes at room temperature.

ABTS Liquid Substrate:

(ABTS Substrate should be at ambient temperature prior to use)

Aspirate and wash plate 4 times. Add 100 μ l of substrate solution to each well. Incubate at room temperature for color development. Monitor color development with an ELISA plate reader at 405 nm with wavelength correction set at 650 nm.

NOTE: Reliable standard curves are obtained when either O.D. readings do not exceed 0.2 units for the zero standard concentrations, or 1.2 units for the highest standard concentration. The plate should be monitored at 5-minute intervals for approximately 50 minutes.

CROSS REACTIVITY

When tested at 50ng/ml the following antigens exhibited 100% cross reactivity:

Rat IFN- γ

When tested at 50ng/ml the following antigens did not exhibit significant cross reactivity:

Human IFN- α , IFN- β , IFN- γ , IFN- λ 1, IFN λ 2, IL-10, IL-18

Murine IFN λ 2, IL-10, IL-12, IL-18

Rat IL-10

