$\mathsf{BD}\;\mathsf{OptEIA}^{\mathsf{TM}}$

Cat. No. 555165

Calculation of Results

Calculate the mean absorbance for each set of duplicate standards, controls and samples. Subtract the mean zero standard absorbance from each.


Plot the standard curve on log-log graph paper, with IL-12 (p40) concentration on the x-axis and absorbance on the y-axis. Draw the best fit curve through the standard points.

To determine the IL-12 (p40) concentration of the unknowns, find the unknown's mean absorbance value on the y-axis and draw a horizontal line to the standard curve. At the point of intersection, draw a vertical line to the x-axis and read the IL-12 (p40) concentration. If samples were diluted, multiply the IL-12 (p40) concentration by the dilution factor.

Computer data reduction may also be employed, utilizing log-log regression analysis.

Typical Standard Curve

This standard curve is for demonstration only. A standard curve must be run with each assav.

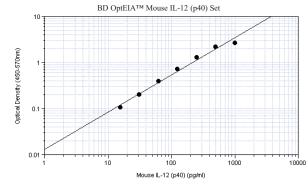

Specificity

Cross Reactivity: The following factors were tested in the BD OptEIATM assay at ≥ 10 ng/mL and no cross-reactivity (value ≥ 4 pg/mL) was identified.

Recombinant Human

IL-1α, IL-1β, IL-2, IL-3, IL-4, IL-5, IL-6, IL-7, IL-8, IL-9, IL-10, IL-11, IL-12 (p40), IL-12 (p70), IL-13, IL-15, G-CSF, GM-CSF, IFN-y, CD23, Lymphotactin, MIP-1α, MIP-1β, MCP-1, MCP-2, NT-3, PDGF-AA, SCF, TNF, LT-α (TNF-β), VEGF

Recombinant Mouse

IL-1β, IL-2, IL-3, IL-4, IL-5, IL-6, IL-7, IL-9, IL-10, IL-15, IFN-γ, GM-CSF, MCP-1, TCA3, TNF

Recombinant Rat IL-2, IL-4, IL-6, IL-10, GM-CSF, IFN-y, TNF

Other:

Viral IL-10 (1 ng/mL), Rabbit TNF

Note: This ELISA set will recognize mouse IL-12 (p70) [i.e it will recognize the (p40) subunit in the context of the (p70) complex]. To account for the complexed subunits, please utilize the BD OptEIATM mouse IL-12 (p70) ELISA Set MN 555256.

Note: Although not directly tested by BD Biosciences, the IL-12 (p40) subunit has been reported to be shared with IL-23.7

Standardization

This immunoassay is calibrated against recombinant mouse IL-12 (p40).

Assay Optimization

- 1. BD OptEIA[™] Sets allow flexible assay design to fit individual laboratory needs. To design an immunoassay with different sensitivity and dynamic range, the following parameters can be varied: Capture, Detection Antibody titers, Incubation time, Incubation temperature, Assay Diluent formulation, Buffer pH, ionic strength, protein concentration, Type of substrate, Washing technique (i.e., number of wash repetitions and soak times)
- 2. "Typical Standard Curve" and 20-plate yield were obtained in the BD Biosciences Pharmingen laboratory, using the recommended procedure and manual plate washing.

Troubleshooting

Poor Precision Possible Source · Inadequate washing/ aspiration of wells · Inadequate mixing of reagents · Imprecise/ inaccurate pipetting · Incomplete sealing of plate

Poor Standard Curve Possible Source · Improper standard handling/ dilution standards · Incomplete washing/ aspiration of wells · Imprecise/ inaccurate pipetting · Improper buffer/ diluent used

Low Absorbances Possible Source · Inadequate reagent volumes added to wells · Incorrect incubation times/ temperature

> · Incorrect antibody titration Detector preparation Improper buffer/ diluent used

· Overly high wash/aspiration pressure from automated plate-washer

Limitations of the Procedure

• Samples that generate absorbance values higher than the standard curve should be diluted with Standard Diluent and re-assayed.

• Interference by drug metabolites, soluble receptors, or other binding proteins in specimens has not been thoroughly investigated. The possibility of interference cannot be excluded.

• BD OptEIATM Sets are intended for use as an integral unit. Do not mix reagents from different Set batches. Reagents from other manufacturers are not recommended for use in this Set.

References

- 1. Chougnet, et al. J Immunol. 2001; 166; 3210-3217.
- 2. Iwasaki A, et al. J Immunol. 2001; 166: 4884-4890.
- 3. Iwasaki A, et al. J Exp Med. 2000; 190: 229-239.
- 4. He J. et al. J Exp Med. 2000; 191: 1605-1610.
- 5. Morelli AE, et al. J Virol. 2000; 74: 9617-9628.
- 6. Ashkar S, et al. Science. 2000; 287: 860-864.
- 7. Oppmann B, et al. Immunity. 2000; 13: 715-725.

Corrective Action · Check function of washing system · Ensure adequate mixing · Check/ calibrate pipettes

· Check function of washing system

Corrective Action · Check/ calibrate pipettes times/reagents warmed to RT

Check buffer/ diluent preparation pH

United States 877.232.8995

Technical Data Sheet

Materials Provided

fied in this package.

Capture Antibody

Detection Antibody

Enzyme Reagent

Standards

(lot-specific)

Anti-Mouse IL-12 monoclonal antibody

Biotinylated Anti-Mouse IL-12 monoclonal antibody

Recombinant mouse IL-12 (p40), lyophilized

Instruction / Analysis Certificate

Streptavidin-horseradish peroxidase conjugate (SAv-HRP)

Mouse IL-12 (p40) ELISA Set

The OptEIA[™] Set for mouse interleukin-12 (p40) (IL-12 (p40)) contains

assays (ELISA) for natural or recombinant mouse IL-12 (p40) in serum,

plasma, and cell culture supernatants. Sufficient materials are provided to

yield approximately 20 plates of 96-wells if the recommended storage,

materials, buffer preparation, and assay procedure are followed as speci-

the components necessary to develop enzyme-linked immunosorbent



Canada 866.979.9408

32.2.400.98.95

0120.8555.90

Asia/Pacific

65.6861.0633

55.11.5185.9995

Latin America/Caribbean

Europe

Japan

BD Biosciences Pharmingen

10975 Torreyana Road San Diego, CA 92121 Customer/Technical Service Tel 877.232.8995 (US) Fax 858.812.8888 www.bdbiosciences.com

Conditions: The information disclosed herein is not to be construed as a recommendation to use the above product in violation of any patents. BD Biosciences will not be held responsible for patent infringement of other violations that may occur with the use of our products. Purchase does not include or carry any right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of Becton Dickinson and Company is strictly prohibited

Unless otherwise noted, BD, BD Logo and all other trademarks are the property of Becton, Dickinson and Company. ©2015 BD

· Ensure complete seal on plate Corrective Action · Ensure correct preparation, storage of

· Check/ calibrate pipettes · Check buffer/ diluent preparation, pH

· Ensure sufficient incubation · Check Capture Ab and Working

· Utilize manual washing

Recommended buffers, solutions

Note: Do not use sodium azide in these preparations. Sodium azide inactivates the horseradish peroxidase enzyme.

The BD OptEIA[™] Reagent Set A (Cat. No 550536) containing Coating Buffer, Assay Diluent, Substrate Reagents A and B, Stop Solution and 20X Wash Buffer Concentrate is recommended.

- Coating Buffer 0.2 M Sodium Phosphate, pH 6.5 12.49 g Na₂HPO₄, 15.47 g NaH₂PO₄; q.s. to 1.0 L; pH to 6.5. Freshly prepare or use within 7 days of preparation, stored at 2-8°C.
- Assay Diluent- PBS* with 10% FBS[#], pH 7.0. The BD Pharmingen[™] Assay Diluent (Cat. No. 555213) is recommended.
 *Phosphate-Buffered Saline: 80.0 g NaCl, 11.6 g Na₂HPO₄, 2.0 g

KH₂PO₄, 2.0 g KCL, q.s. to 10 L; pH to 7.0.

#Fetal Bovine Serum: Hyclone Cat. No. SH30088 (heatinactivated) recommended. Freshly prepare or use within 3 days of preparation, with 2-8°C storage.

 Wash Buffer - PBS* with 0.05% Tween-20. Freshly prepare or use within 3 days of preparation, stored at 2-8°C.

 Substrate Solution - Tetramethylbenzidine (TMB) and Hydrogen Peroxide. The BD Pharmingen[™] TMB Substrate Reagent Set (Cat. No. 555214) is recommended.

5. Stop Solution - 1 M H₃PO₄ or 2 N H₂SO₄

Additional Materials Required

- 96-well Nunc-ImmunoTM polystyrene Maxisorp ELISA flat bottom plates (ThermoFisher Scientific Cat. No. 442404) are recommended
- 2. Microplate reader capable of measuring absorbance at 450 nm
- 3. Precision pipettes
- 4. Graduated cylinder, one liter
- 5. Deionized or distilled water
- 6. Wash bottle or automated washer
- 7. Log-log graph paper or automated data reduction
- 8. Tubes to prepare standard dilutions
- 9. Laboratory timer
- 10. Plate sealers or parafilm

Storage Information

- 1. Store unopened reagents at 2-8°C. Do not use reagents after expiration date, or if turbidity is evident.
- 2. Before use, bring all reagents to room temperature (18-25°C). Immediately after use, return to proper storage conditions.
- 3. Lyophilized standards are stable until expiration date. See below for reconstituted standard storage information.

Specimen Collection and Handling

Specimens should be clear, non-hemolyzed and non-lipemic.

Cell culture supernatants: Remove any particulate material by centrifugation and assay immediately or store samples at \leq -20°C. Avoid repeated freeze-thaw cycles.

Serum: Use a serum separator tube and allow samples to clot for 30 minutes, then centrifuge for 10 minutes at 1000 x g. Remove serum and assay immediately or store samples at \leq -20° C. Avoid repeated freeze-thaw cycles.

Plasma: Collect plasma using citrate, EDTA, or heparin as anticoagulant. Centrifuge for 10 minutes at 1000 x g within 30 minutes of collection. Assay immediately or store samples at \leq -20° C. Avoid repeated freeze-thaw cycles.

Standards Preparation and Handling


- Reconstitution: After warming lyophilized standard to room temperature, carefully open vial to avoid loss of material. Reconstitute lyophilized standard with 1.0 mL of deionized water to yield a stock standard. Allow the standard to equilibrate for at least 15 minutes before making dilutions. Vortex gently to mix.
- Storage/ handling of reconstituted standard: After reconstitution, immediately aliquot standard stock in polypropylene vials at 50 µl per vial and freeze at -80°C for up to 6 months. If necessary, store at 2-8° C for up to 8 hours prior to aliquotting/freezing. Do not leave reconstituted standard at room temperature.

3. Standards Preparation for Assay:

a. Prepare a 1000 pg/mL standard from the stock standard. Vortex to mix. (See dilution instructions on Instruction/Analysis Certificate.)
b. Add 300 μL Assay Diluent to 6 tubes. Label as 500 pg/mL, 250 pg/mL, 125

b. Add 300 µL Assay Diluent to 6 tubes. Label as 500 pg/mL, 250 pg/mL, 1 pg/mL, 62.5 pg/mL, 31.3 pg/mL, and 15.6 pg/mL.

c. Perform serial dilutions by adding 300 μ L of each standard to the next tube and vortexing between each transfer. Assay Diluent serves as the zero standard (0 pg/mL).

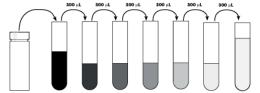

Stoek Standard 1000 pg/ml. 500 pg/mL 250 pg/mL 125 pg/mL 62.5 pg/mL 31.3 pg/mL 15.6 pg/mL

Serial dilutions within the plate may also be performed by pipetting 100 μL of Assay Diluent into each standard well except the highest (1000 pg/mL), then adding 100 μL of the 1000 pg/mL standard to both that well and the 500 pg/mL well, mixing the well contents by rinsing the pipette tip, and adding 100 μL of the 500 pg/mL standard to the 250 pg/mL vell. Continue these dilutions to the 15.6 pg/mL standard well, out of which the extra 100 μL should be discarded.

Working Detector Preparation

(Note: One-step incubation of Biotin/Streptavidin reagents.) Add required volume of Detection Antibody to Assay Diluent. Within 15 minutes prior to use, add required quantity of Enzyme Reagent, vortex or mix well. For recommended dilutions, see lot-specific Instruction/Analysis Certificate. For a full 96-well plate, prepare 12 mL of Working Detector. Discard any remaining Working Detector after use.

Warnings and Precautions

- 1. Reagents which contain preservatives may be toxic if ingested, inhaled, or in contact with skin.
- 2. Handle all serum and plasma specimens in accordance with NCCLS guidelines for preventing transmission of blood-borne infections.
- Capture Antibody contains < 0.1% sodium azide. Sodium azide yields highly toxic hydrazoic acid under acidic conditions. Dilute azide compounds in running water before discarding to avoid accumulation of potentially explosive deposits in plumbing.
- 4. Detection Antibody contains BSA and ProClin® -150 as a preservative.
- 5. Enzyme Reagent contains BSA and ProClin® -150 as a preservative.
- 6. Source of all serum proteins is from USDA inspected abattoirs located in the United States.
- Warning: Recombinant Mouse IL-12 (p40) lyophilized standard (component 51-26196E) and Detection Antibody Biotin Anti-Mouse IL-12 (p40) (component 51-26192E) contains 0.02% (w/w) of a CMIT/MIT mixture (3:1), which is a mixture of: 5-chloro-2-methyl-4-isothiazolin-3-one [EC No 247-500-7] and 2methyl-4-isothiazolin-3-one [EC No 220-239-6] (3:1).

Hazard statements

May cause an allergic skin reaction. *Precautionary statements*

- Wear protective gloves / eye protection.
- Wear protective clothing.
- Avoid breathing mist/vapours/spray.

If skin irritation or rash occurs: Get medical advice/attention.

IF ON SKIN: Wash with plenty of water.

Dispose of contents/container in accordance with local/regional/national/interna-

tional regulations.

Recommended Assay Procedure

- Coat microwells with 100 μL per well of Capture Antibody diluted in Coating Buffer. For recommended antibody coating dilution, see lot-specific Instruction/Analysis Certificate. Seal plate and incubate overnight at 4° C.
- Aspirate wells and wash 3 times with ≥ 300 μL/well Wash Buffer. After last wash, invert plate and blot on absorbent paper to remove any residual buffer.
- 3. Block plates with $\geq 200 \ \mu$ L/well Assay Diluent. Incubate at RT for 1 hour.
- 4. Aspirate/wash as in step 2.
- 5. Prepare standard and sample dilutions in Assay Diluent. See "Standards Preparation and Handling".
- Pipette 100 μL of each standard, sample, and control into appropriate wells. Seal plate and incubate for 2 hours at RT.
- 7. Aspirate/ wash as in step 2, but with 5 total washes.
- Add 100 µL of Working Detector (Detection Antibody + SAv-HRP reagent) to each well. Seal plate and incubate for 1 hour at RT.
- 9. Aspirate/ wash as in step 2, but with 7 total washes. Note: In this final wash step, soak wells in wash buffer for 30 seconds to 1 minute for each wash.
- 10. Add 100 μ L of Substrate Solution to each well. Incubate plate (without plate sealer) for 30 minutes at room temperature in the dark.
- 11. Add 50 μL of Stop Solution to each well.
- Read absorbance at 450 nm within 30 minutes of stopping reaction. If wavelength correction is available, subtract absorbance at 570 nm from absorbance 450 nm.

Assay Procedure Summary

- 1. Add 100 µL diluted Capture Ab to each well.
 - Incubate overnight at 4°C.
- 2. Aspirate and wash 3 times.
- 3. Block plates: 200 µL Assay Diluent to each well. Incubate 1 hr RT
- 4. Aspirate and wash 3 times.
- 5. Add 100 μ L standard or sample to each well. Incubate 2 hr RT.
- 6. Aspirate and wash 5 times.
- Add 100 μL Working Detector (Detection Ab + SAv-HRP) to each well. Incubate 1 hr RT
- 8. Aspirate and wash 7 times (with 30 sec. to 1 min soaks)
- 9. Add 100 μ L Substrate Solution to each well. Incubate 30 min RT in dark
- 10. Add 50 μL Stop Solution to each well. Read at 450 nm within 30 min with λ correction 570 nm.