

UltraPure™ Glycogen

Cat. no. 10814-010

Size 100 µL Store at -30°C to -10°C

Doc. Part no. 10814010.pps

Pub. no. MAN0000908

Rev. 2.0

Description

Use UltraPure™ glycogen as a carrier for nucleic acid precipitation in ethanol at a concentration of $0.2-0.4 \,\mu\text{g}/\mu\text{L}$ of aqueous sample. UltraPure $^{\text{TM}}$ glycogen is supplied at 20 µg/µL in RNase-free water.

Notes

- Use UltraPure[™] glycogen as a carrier in ethanol precipitation of ligation reactions. Using 8 µg of UltraPure™ glycogen in a 100-µL reaction with chemically competent cells does not affect the transformation (in-house communication).
- Used at a concentration of 0.001 mg/mL, UltraPure[™] glycogen does not inhibit TdT.
- A concentration of ≤ 2 mg/mL glycogen has no effect on SuperScript® II Reverse Transcriptase (Gerard, 1994).
- T4 RNA Ligase is not inhibited by samples precipitated in the presence of 0.02 mg/mL glycogen (Apte, 1993).
- UDG is not inhibited by glycogen.
- Glycogen can interfere with DNA-protein interactions (Galliard, 1990).
- Glycogen can decrease RNA quality and inhibit RT-PCR in RNA isolated from tissues rich in polysaccharides and proteoglycans (Chomczynski, 1995).

For Research Use Only. Not for use in diagnostic procedures.

References

Apte, A. N. and Siebert, P. D. (1993) Biotechniques 15.5, 890.

Chomcyznski, P. and Mackey, K. (1995) Biotechniques 19.6, 942.

Galliard, C. and Strauss, F. (1990) Nucleic Acids Research 18, 378.

Gerard, G. (1994) Focus® 16:4, 102.

Certificate of Analysis

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to **www.lifetechnologies.com/support** and search for the Certificate of Analysis by product lot number, which is printed on the box.

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

©2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

DISCLAIMER - LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT. INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

For support visit lifetechnologies.com/support or email techsupport@lifetech.com

lifetechnologies.com

life technologies™