

SuperScript[®] VILO[™] MasterMix

Part no. 100012386

MAN0004286

Rev. Date: 6 June 2011

Cat. nos:

11755-050

11755-250

11755-500

Size:

50 reactions

250 reactions

500 reactions

Store at -20°C (non-frost-free)

Description

SuperScript[®] VILO[™] MasterMix provides the high-temperature capability of SuperScript[®] III Reverse Transcriptase (RT) in an optimized format for generating first-strand cDNA for use in real-time quantitative RT-PCR (qRT-PCR). This formulation can be used with very low and very high amounts of input RNA (up to 2.5 µg total RNA in a 20-µL reaction).

SuperScript[®] VILO[™] Master Mix includes SuperScript[®] III RT, RNaseOUT[™] Recombinant Ribonuclease Inhibitor, a proprietary helper protein, random primers, MgCl₂, and dNTPs.

SuperScript[®] III RT is an engineered version of M-MLV RT with reduced RNase H activity and increased thermal stability. The enzyme can be used to synthesize cDNA at a temperature range of 42–55°C. Because SuperScript[®] III RT is not significantly inhibited by ribosomal and transfer RNA, it can be used to synthesize cDNA from total RNA. RNaseOUT[™] Recombinant Ribonuclease Inhibitor safeguards against the degradation of target RNA due to ribonuclease contamination.

Component	50 rxn kit	250 rxn kit	500 rxn kit
SuperScript [®] VILO [™] MasterMix	200 µL	1000 µL	2000 µL

Intended Use: For research use only.

Not intended for any animal or human therapeutic or diagnostic use.

Guidelines for cDNA Synthesis

- High-quality, intact RNA is essential for accurate quantification in qPCR. RNA should be devoid of RNase contamination and aseptic conditions should be maintained. RNA quality can be analyzed using a bioanalyzer or by agarose gel electrophoresis.
- Starting material can range up to 2.5 µg of total RNA in a 20-µL cDNA synthesis reaction. RNA quantity can be determined using UV absorbance at 260 nm or the Qubit® RNA Assay Kit and Qubit® 2.0 Fluorometer (see page 4 for ordering information).
- To isolate total RNA, we recommend TRIzol® Reagent, the PureLink® RNA Mini Kit, or the MagMAX™-96 Total RNA Isolation Kit (see page 4). Isolation of mRNA is typically not necessary, although incorporating this step may improve the yield of specific cDNAs.
- DNase I, Amplification Grade, may be used to eliminate genomic DNA contamination from the total RNA (see page 4).
- Shorter incubation times and/or higher temperatures may be used (e.g., 50°C for 30 minutes), but may result in reduced yields of cDNA.
- For increased yields of cDNA, longer incubation times may be used (up to 120 minutes at 42°C).

qPCR Using Fluorescent Primers or Probes

Up to 10% of the qPCR reaction volume may be undiluted cDNA (e.g., for a 20-µL qPCR, use up to 2 µL of undiluted cDNA).

qPCR Using SYBR® Green or SYBR® GreenER™ Reagent

If you started with ≤100 ng of total RNA, up to 10% of the qPCR reaction volume may be undiluted cDNA (e.g., for a 20-µL qPCR, use up to 2 µL of undiluted cDNA).

If you started with >100 ng total RNA, we recommend testing a serial dilution of cDNA in qPCR for optimal results. Higher concentrations of cDNA may affect the signal baseline in SYBR® Green and SYBR GreenER™ reactions.

First-Strand cDNA Synthesis

The following protocol has been optimized for generating first-strand cDNA for use in two-step qRT-PCR. The reaction volume may be scaled as needed up to 100 μL . A negative RT control protocol is provided below.

1. For a single reaction, combine the following components in a sterile PCR tube or plate well on ice.

Component	Volume
SuperScript [®] VILO [™] MasterMix	4 μL
RNA (up to 2.5 μg)	X μL
DEPC-treated water	to 20 μL

2. Gently mix and incubate at 25°C for 10 minutes.
3. Incubate at 42°C for 60 minutes.
4. Terminate the reaction at 85°C at 5 minutes.
5. Use the diluted or undiluted cDNA in qPCR or store at -20°C.

Negative RT Control

1. For a volume of RNA = X μL , add the following to a sterile PCR tube or plate well on ice.

Component	Volume
SuperScript [®] VILO [™] MasterMix	4 μL
DEPC-treated water	16 - X μL

2. Incubate at 65°C for 10 minutes to denature the reverse transcriptase.
3. Add X μL of RNA (up to 2.5 μg) for a total reaction volume of 20 μL .
4. Proceed with steps 2-5 from the "First-Strand cDNA Synthesis" protocol above.

Product Qualification

The Certificate of Analysis (CofA) provides detailed quality control information for each product. The CofA is available on our website at www.invitrogen.com/cofa, and is searchable by product lot number, which is printed on each box.

Additional Products

	Amount	Catalog number
TRIzol® Reagent	100 ml	15596-026
	200 ml	15596-018
PureLink® RNA Mini Kit	10 preps	12183020
	50 preps	12183018A
MagMAX™-96 Total RNA Isolation Kit	96 rxns	AM1830
Qubit® RNA Assay Kit	500 assays	Q32855
Qubit® 2.0 Fluorometer	1 unit	Q32866
DNase I, Amplification Grade	100 units	18068-015

Limited Use Label License No. 358: Research Use Only

The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

©2011 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners. TRIzol is a registered trademark of Molecular Research Center, Inc.

For support visit www.lifetechnologies.com/support or email techsupport@lifetech.com

www.lifetechnologies.com

