

Platinum[®] PCR SuperMix High Fidelity

Cat. No. 12532-016

Size: 100 Reactions

Cat. No. 12532-024

Size: 5000 Reactions

Store at -20°C in a
non-frost-free freezer.

Description:

Platinum[®] PCR SuperMix High Fidelity provides qualified reagents for high fidelity amplification of DNA templates by polymerase chain reaction (PCR). It is effective over a large range of target sizes—up to 15 kb genomic DNA. The mixture contains anti-*Taq* DNA polymerase antibody, Mg⁺⁺, deoxyribonucleotide triphosphates, recombinant *Taq* DNA polymerase, and *Pyrococcus* species *GB-D* thermostable polymerase. Anti-*Taq* DNA polymerase antibody inhibits polymerase activity, providing an automatic “hot start” (1,2) and permitting room temperature set-up (polymerase activity is restored after a denaturation step in PCR cycling at 94°C). Antibody-mediated hot starts improve PCR specificity and yield (3). *Pyrococcus* species *GB-D* polymerase possesses a proofreading ability by virtue of its 3' to 5' exonuclease activity (3). Mixture of the proofreading enzyme with *Taq* DNA polymerase increases fidelity approximately six times over that of *Taq* DNA polymerase alone.

Platinum[®] PCR SuperMix High Fidelity is supplied at 1.1X concentration to allow approximately 10% of the final reaction volume to be used for the addition of primer and template solutions. Reagents sufficient for 100 or 5000 amplification reactions of 50 µl each are provided.

Component:

Platinum[®] PCR SuperMix High Fidelity*

100 Rxn Kit

4 × 1.125 ml

5000 Rxn Kit

4 × 56.25 ml

*22 U/ml complexed recombinant *Taq* DNA polymerase, *Pyrococcus* species *GB-D* thermostable polymerase, and Platinum[®] *Taq* Antibody; 66 mM Tris-SO₄ (pH 8.9); 19.8 mM (NH₄)₂SO₄; 2.4 mM MgSO₄; 220 µM dNTPs; and stabilizers.

Part. No. 12532.pps

Rev. date: 11 May 2010

Storage Conditions:

After thawing, Platinum® PCR SuperMix High Fidelity may be stored at 4°C for 3 months or -20°C for 1 year. Storage at 4°C avoids the necessity of thawing the mix before assembling the reaction. There is no detectable decrease in enzyme activity or performance after storage for 3 months at 4°C, or after 15 freeze-thaw cycles.

Recommendations and Guidelines:

- Because PCR is a powerful technique capable of amplifying trace amounts of DNA, take all appropriate precautions to avoid cross-contamination. Ideally, amplification reactions should be assembled in a DNA-free environment.
- Reactions may be assembled either at room temperature or on ice. We have observed no significant difference in reaction efficiency between these setup conditions.
- For multiple reactions, you can prepare a master mix of Platinum® PCR SuperMix High Fidelity and the component(s) common to all reactions.

PCR Protocol:

1. Add the following components in any order to each reaction tube:
 - a. 45 µl Platinum® PCR SuperMix High Fidelity
 - b. Primer solution (200 nM final concentration of each is recommended)*
 - c. Template DNA solution (1–200 ng genomic DNA)*

*A standard 50-µl PCR reaction includes a combined primer and template volume of 5 µl; we have observed no decrease in product yield if the amount of primer and template solution is between 1 µl and 15 µl.
2. Mix contents of tubes and cover with mineral or silicone oil if necessary.
3. Cap tubes and load in thermal cycler.
4. Incubate tubes at 94°C for 30 s to 2 min to completely denature the template and activate the enzyme.
5. Perform 25-35 cycles of PCR amplification as follows:

Denature	94°C for 15-30 s
Anneal	55°C for 15-30 s
Extend	68°C for 1 min per kb

Tips:

- If the PCR efficiency is not optimal, repeat the reaction with different primer concentrations from 100 to 500 nM, in 100 nM increments.
- For longer genomic DNA targets (>15 kb), we recommend adding 1–1.5 U of Platinum® *Taq* DNA Polymerase (Cat. no. 10966-018) to the reaction mix and increasing the extension time as specified (1 min per kb).
- At higher volumes of primer and template, the MgSO₄ concentration in the reaction will drop to suboptimal levels and yield will decrease. For combined primer-template volumes of >15 µl (in solution with 45 µl of Platinum® PCR SuperMix High Fidelity), we recommend adjusting the final MgSO₄ concentration in the reaction to 2.2 mM.

Quality Control:

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to www.invitrogen.com/support and search for the Certificate of Analysis by product lot number, which is printed on the box.

References:

1. Chou, Q., Russel, M., Birch, D., Raymond, J., Bloch, W. (1992) *Nucl. Acids Res.*, 20, 1717.
2. Sharkey, D.J., Scalice, E.R., Christy, K.G., Atwood, S.M., Daiss, J.L. (1994) *BioTechnology*, 12, 506.
3. Westfall, B., Sitaraman, K., Solus, J., Hughes, J., Rashtchian, A. (1997) *Focus*®, 19,2, 46.

Limited Use Label License No. 1: Thermostable Polymerases

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,789,224, 5,618,711, and 6,127,155. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim, no right to perform any patented method, and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

Limited Use Label License No: 5 Invitrogen Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. For products that are subject to multiple limited use label licenses, the terms of the most restrictive limited use label license shall control. Life Technologies Corporation will not assert a claim against the buyer of infringement of patents owned or controlled by Life Technologies Corporation which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Life Technologies is willing to accept return of the product with a full refund. For information about purchasing a license to use this product or the technology embedded in it for any use other than for research use please contact Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008 ; Phone (760) 603-7200 or e-mail: outlicensing@lifetech.com

Limited Use Label License No: 14 Direct Inhibition by Anti-polymerase Antibodies

Licensed to Life Technologies Corporation, under U.S. Patent Nos. 5,338,671; 5,587,287, and foreign equivalents for use in research only.

Limited Use Label License No: 117 Thermostable DNA Polymerase Blend

This product is the subject of U.S. Patent No. 5,436,149 and corresponding foreign patents licensed by Takara Corporation to Life Technologies Corporation. This product may be used only for research purposes and may not be resold or used for therapeutic purposes.

©2010 Life Technologies Corporation. All rights reserved.

For research use only. Not intended for any animal or human or diagnostic use.