

# Human Mammary Epithelial Cells (HMEC)

Cat. no. A10565

Quantity:  $\geq 500,000$  viable cells/vial

## Product Description

HMEC are normal human mammary epithelial cells isolated from reduction mammaplasty tissue. Each vial of this product contains  $\geq 5 \times 10^5$  viable cells are cryopreserved at the end of the 6' culture in a medium containing 10% DMSO. Approved lots of cells have tested positive by immunofluorescent methods for cytokeratins 5/6, 8, 18, and E-cadherin. An independent laboratory tests the cells for the presence of mycoplasma, Hepatitis B, Hepatitis C, and HIV-1 viruses. These agents were not detected. Each lot of cells is performance tested by culturing the cells through multiple passages in complete HuMEC medium in the absence of antibiotics and antimycotics. During this culture period, no contamination by bacteria, yeast, or other fungi was detected. Upon thawing, the cells are guaranteed to be  $\geq 70\%$  viable (by trypan blue staining) and have a potential of  $\geq 16$  population doublings when handled according to the directions provided in this document. For recommended precautions to be taken when handling human cells, please read the caution statement.

## Intended Use

Cryopreserved HMEC are intended for use by researchers investigating the molecular and biochemical basis of various normal and disease processes. **This product is for research use only. Not intended for human or animal therapeutic or diagnostic use.**

## Storage and Stability

Cryopreserved HMEC should arrive frozen on dry ice. If the cells are not to be used immediately, prepare a space for storage of the vial in the vapor phase of a liquid nitrogen freezer. While wearing protective eyewear, gloves, and a laboratory coat, remove the vial from its shipping container and place immediately in the liquid nitrogen freezer. Although the viability of cryopreserved cells decreases with time in storage, useful cultures can usually be established even after 2 years of storage at liquid nitrogen temperatures.

## Caution

Although cryopreserved cells have been tested for the presence of various hazardous agents, diagnostic tests are not necessarily 100% accurate. In addition, human cells may harbor other known or unknown agents, or organisms which could be harmful to your health or cause fatal illness. Treat all human cells as potential pathogens. Wear protective clothing and eyewear. Practice appropriate disposal techniques for potentially pathogenic or biohazardous materials. In case of contact with eyes, rinse immediately with plenty of water and seek medical advice.

## Initiating Cultures from Cryopreserved Cells

HMEC Culture System from Invitrogen		
Description	Cat. no.	Size
HuMEC Ready Medium Contains:	12752-010	1 kit
HuMEC Basal Serum Free Medium	12753-018	1 x 500 ml
HuMEC Supplement	12754-016	1 x 5 ml
Bovine Pituitary Extract	13028-014	1 x 25 mg
Trypsin/EDTA	R-001-100	100 ml
Trypsin Neutralizer	R-002-100	100 ml

We recommend seeding HMEC recovered from cryopreservation at a density of  $2.5 \times 10^3$  viable cells/cm<sup>2</sup>. For example,  $3 \times 75$  cm<sup>2</sup> or  $9 \times 25$  cm<sup>2</sup> tissue culture flasks are usually established from one vial containing  $\geq 5 \times 10^5$  HMEC. The procedure given below is a sample protocol for establishing cultures from one vial.

1. Prepare a bottle of supplemented HuMEC medium according to the instructions supplied with that product.
  2. Remove a vial of HMEC from liquid nitrogen storage, taking care to protect hands and eyes.
  3. Dip the lower half of the vial into a 37°C water bath to thaw.
  4. When the contents of the vial have just thawed, wipe the outside of the vial with disinfecting solution and move to a Class II, type A laminar flow culture hood.
  5. Open the vial and gently pipet the suspension up and down with a 1 ml pipette to disperse the cells.
  6. Remove 20  $\mu$ l from the vial and dilute the cell suspension in 20  $\mu$ l of trypan blue solution (for example: Invitrogen Trypan Blue Stain Cat. no. 15250-061).
  7. Use a hemacytometer to determine the number of viable cells per ml.
- Note:** Do not warm the reagents prior to use.
8. Dilute the contents of the vial (1 ml) to a concentration of  $1.25 \times 10^4$  viable cells/ml using the supplemented medium from step 1, above.
  9. Add 5 ml of cell suspension to each 25 cm<sup>2</sup> culture flask or 15 ml of cell suspension to each 75 cm<sup>2</sup> culture flask.
  10. Following inoculation, swirl the medium in the flasks to evenly distribute the cells.
  11. Incubate the cultures in a 37°C, 5% CO<sub>2</sub>/95% air, humidified cell culture incubator. For best results, do not disturb the culture for at least 18 hours after the culture has been initiated.

**For research use only.**

Life Technologies Corporation • 5791 Van Allen Way • Carlsbad • CA 92008 • Tel: 800.955.6288 • [www.invitrogen.com](http://www.invitrogen.com) • E-mail: tech\_support@invitrogen.com

## Maintaining Stock Cultures

1. Change the culture medium to freshly supplemented medium, 18 to 24 hours following plating of cryopreserved cells. For subsequent subcultures, change the medium 48 hours after establishing the subculture.
2. Change the medium every other day thereafter, until the culture is approximately 50% confluent.
3. Once the culture reaches 50% confluency, change the medium every day until the culture is approximately 80-90% confluent.

### Notes

- To achieve the highest cell densities, change the culture medium every day as the cultures approach confluency. To obtain rapidly proliferating subcultures, subculture HMEC before they become more than 90% confluent. The number of subcultures (passages) that can be achieved vary with the starting cell density and the methods employed by individual investigators.
- HMEC cultures seeded at  $2.5 \times 10^3$  cells/cm<sup>2</sup> from cryopreserved cells should reach 80-90% confluency in 5-7 days. At this time, most of the cells should have epithelial or "cobblestone" morphology. Some irregularly sized and shaped cells may be observed.

## Subculture of HMEC

View the culture under the microscope to confirm that it is subconfluent (80-90%), and that there are mitotic cells present. This protocol is designed for the subculture of one 25 cm<sup>2</sup> culture flask. If different-sized culture vessels are used, adjust the reagent volumes accordingly.

1. Assemble supplemented medium and subculture reagents according to the table on page 1.

**Note:** Do not warm the reagents prior to use.

2. Assemble the appropriate culture vessels, sterile pipettes, and sterile 15 ml conical tubes.
3. Remove all of the culture medium from the flask.
4. Add 3 ml of Trypsin/EDTA solution to the flask. Rock the flask to ensure that the entire surface is covered.
5. Immediately remove all 3 ml of Trypsin/EDTA solution from the flask.

### Limited Use Label License No. 5: Invitrogen Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. For products that are subject to multiple limited use label licenses, the terms of the most restrictive limited use label license shall control. Life Technologies Corporation will not assert a claim against the buyer of infringement of patents owned or controlled by Life Technologies Corporation which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Life Technologies is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Life Technologies Corporation, 5791 Van Allen Way, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500. Email: [outlicensing@invitrogen.com](mailto:outlicensing@invitrogen.com). ©2009 Life Technologies Corporation. All rights reserved.

**For research use only.**

Life Technologies Corporation • 5791 Van Allen Way • Carlsbad • CA 92008 • Tel: 800.955.6288 • [www.invitrogen.com](http://www.invitrogen.com) • E-mail: [tech\\_support@invitrogen.com](mailto:tech_support@invitrogen.com)

6. Add 1 ml of fresh Trypsin/EDTA solution to the flask.
7. Incubate the flask at 37°C for 4-6 minutes. View the culture under a microscope to confirm cells have become round.
8. Rap the flask very gently to dislodge cells from the surface of the flask.
9. Add 3 ml of Trypsin Neutralizer solution to the flask and transfer the detached cells to a sterile 15 ml tube.
10. Add 3 ml of additional Trypsin Neutralizer solution to the flask and pipette the solution over the flask surfaces several times to remove any remaining cells. Add this suspension to the 15 ml conical tube.
11. Centrifuge the cells at 180 × g for 7 minutes. Observe the cell pellet.
12. Remove the supernatant from the tube, being careful not to dislodge the cell pellet.
13. Resuspend the cell pellet in 4 ml supplemented medium. Pipet the cells up and down with a 10 ml pipette to ensure a homogeneous cell suspension.
14. Use a hemacytometer to determine the number of viable cells per ml.
15. Dilute the cells in supplemented medium and seed new culture vessels at  $2.5 \times 10^3$  viable cells/cm<sup>2</sup>.
16. Incubate the cultures in a 37°C, 5% CO<sub>2</sub>/95% air, humidified cell culture incubator.

### Note

Damage to cultured HMEC can occur during trypsinization. This damage may result from exposure of the cells to the Trypsin/EDTA solution for excessive lengths of time and/or excessive mechanical agitation. HMEC cultures can alternatively be trypsinized at room temperature incubation for 10-12 minutes. Monitor cell detachment and otherwise follow "Subculture of HMEC" above. Invitrogen TrypLE™ Express (Cat. no. 12604-013) may be substituted for Trypsin/EDTA without affecting HMEC performance. Invitrogen Penicillin/Streptomycin (Cat. no. 15140-148), and Gentamicin/Amphotericin (Cat. no. R-015-10) reagents have been performance validated for routine use and subculture with HMEC.