

# FreeStyle™ 293-F Cells

**Catalog Number** R790-07

**Document Part Number** 25-0457

**Publication Number** MAN0000255

**Revision** A.0

Information in this document is subject to change without notice.

## **DISCLAIMER**

TO THE EXTENT ALLOWED BY LAW, LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

## **LIMITED USE LABEL LICENSE No. 561: Internal Research and Development Use of Mammalian and Insect Cell Lines**

Notice to Purchaser: The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product and progeny, to perform internal research and development for the sole benefit of the purchaser. The purchase of this product does not grant the purchaser any additional rights, including, without limitation: (a) the right to use the product or its components in commercial applications of any kind, including bioproduction (e.g. use of the product to manufacture therapeutic agents or diagnostic test components, such as proteins, antibodies, viral particles or virus-like particles), quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration; (b) the right to use the product or its components as a therapeutic agent or diagnostic test component; (c) the right to use the product or its components to produce material for use in human clinical trials; or (d) the right to transfer or resell the product or its components in any form, progeny or derivative. For information on obtaining additional rights, please contact [outlicensing@lifetech.com](mailto:outlicensing@lifetech.com) or Out Licensing, Life Technologies Corporation, 5791 Van Allen Way, Carlsbad, California 92008.

© 2015 Thermo Fisher Scientific Inc. All rights reserved. Pluronic is a registered trademark of BASF Corporation. Celligen is a registered trademark of New Brunswick Scientific, Co., Inc. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries.

# Contents

|                                |           |
|--------------------------------|-----------|
| Contents and storage.....      | 4         |
| <b>Introduction .....</b>      | <b>5</b>  |
| Overview .....                 | 5         |
| <b>Methods .....</b>           | <b>8</b>  |
| Important guidelines.....      | 8         |
| Thaw and establish cells ..... | 10        |
| Subculture cells .....         | 12        |
| Cryopreservation.....          | 14        |
| Transfect cells .....          | 16        |
| Troubleshooting.....           | 20        |
| <b>Appendix .....</b>          | <b>22</b> |
| Accessory products .....       | 22        |
| Technical support .....        | 23        |
| Purchaser notification .....   | 24        |
| References.....                | 25        |

# Contents and storage

---

## Shipping and storage

This manual is shipped with FreeStyle™ 293-F Cells. FreeStyle™ 293-F Cells are shipped on dry ice. Upon receipt, store in **liquid nitrogen**.

---

## Contents

### **Storage conditions: Liquid nitrogen**

Amount supplied: One vial containing  $1 \times 10^7$  cells

Composition: 1 mL of cells in 90% FreeStyle™ 293 Expression Medium and 10% DMSO.

---



**CAUTION!** Handle as potentially biohazardous material under at least Biosafety Level 2 containment.

**WARNING! GENERAL CHEMICAL HANDLING.** For every chemical, read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support). This product contains Dimethyl Sulfoxide (DMSO); components of the product may be absorbed into the body through the skin.

---

# Introduction

## Overview

---

### Introduction

The FreeStyle™ 293-F cell line is derived from the 293 cell line (see **Parental Cell Line**) and is intended for use with the FreeStyle™ MAX 293 Expression System (Cat. no. K9000-10) or the FreeStyle™ 293 Expression System (Cat. no. K9000-01). FreeStyle™ 293-F cells are adapted to suspension culture in FreeStyle™ 293 Expression Medium. Frozen cells are supplied in and may be thawed directly into FreeStyle™ 293 Expression Medium (see **Thaw and establish cells**, page 10).

---

### Parental cell line

The 293 cell line is a permanent line established from primary embryonal human kidney transformed with sheared human adenovirus type 5 DNA (Graham et al., 1977; Harrison et al., 1977). The E1A adenovirus gene expressed in these cells participates in transactivation of some viral promoters, allowing these cells to produce very high levels of protein.

The FreeStyle™ 293-F cell line is a variant of the 293 cell line that has been adapted to suspension growth in FreeStyle™ 293 Expression Medium. The 293-F cell line was obtained from Robert Horlick at Pharmacopecia.

---

### FreeStyle™ 293-F Cells

Prepared from low passage Master Cell Bank cultures derived from parental 293-F cells that were re-cloned by limiting dilution. The 293 clonal derived cultures are maintained in serum-free conditions for only 30–35 total passages.

- 293-F Master Cell Bank cultures have been tested by an independent service and found to be negative for HBV, HCV, HTLV-I & -II, and HIV-1 & -2
  - Adapted to high density, serum-free, suspension growth and can be maintained in FreeStyle™ 293 Expression Medium.
  - Demonstrates high transfection efficiencies with FreeStyle™ MAX reagent.
  - Suspension cultures may be transfected in FreeStyle™ 293 Expression Medium without the need to change media.
  - Permits transfection of cells at large volumes.
-

## FreeStyle™ 293 Expression Medium

We recommend using FreeStyle™ 293 Expression Medium to grow, maintain, and transfect FreeStyle™ 293-F cells. FreeStyle™ 293 Expression Medium available from Thermo Fisher Scientific (see page 22 for ordering information) is a defined, serum-free formulation specifically developed for the high density, suspension culture and transfection of 293 cells. The medium contains no human or animal origin components and is formulated with Glutamax™-I to increase stability and maximize shelf life. For more information, see our website ([www.lifetechnologies.com](http://www.lifetechnologies.com)) or call Technical Support (see page 23).

---

## Growth of FreeStyle™ 293-F Cells in the medium

Typically, FreeStyle™ 293-F cells cultured in FreeStyle™ 293 Expression Medium demonstrate the following:

- Doubling time in the range of 22–28 hours (doubling time can exceed 28 hours during the first few passages after the cells have been thawed. However, cells may perform poorly if the doubling time is consistently greater than 30 hours.)
- Cell densities of up to  $3 \times 10^6$  cells/mL in shaker or spinner culture
- Cell densities of up to  $4 \times 10^6$  cells/mL in bioreactor culture

Do not allow FreeStyle™ 293-F cell cultures to approach the plateau of culture density before transfection, as this will result in a decrease of transfection efficiency and protein expression.

**Note:** Individual culturing and passaging techniques coupled with cellular heterogeneity inherent within the FreeStyle™ 293-F cell population may result in experimental variability.

---

## FreeStyle™ MAX Reagent

FreeStyle™ MAX Reagent is a proprietary, animal origin-free formulation for the highly efficient transfection of plasmid DNA into eukaryotic cells. The main benefits of FreeStyle™ MAX Reagent are:

- Specifically formulated to achieve the **highest expression levels** and **lowest cytotoxicity** in suspension FreeStyle™ 293-F Cells and FreeStyle™ CHO-S Cells
- No media change required
- Complex formation can be performed in animal-origin free OptiPro™ SFM

For more information, visit [www.lifetechnologies.com](http://www.lifetechnologies.com) or call Technical Support (see page 23).

---

## FreeStyle™ MAX 293 Expression System

The FreeStyle™ 293-F Cells are part of the FreeStyle™ MAX 293 Expression System. Using the FreeStyle™ MAX 293 Expression System for protein production in mammalian cells provides the following advantages:

- Uses 293 human embryonal kidney cells, for very high expression levels of recombinant protein in human cells
- The FreeStyle™ MAX Reagent offers high recombinant protein yield with low cytotoxicity
- Provides a rapid transient transfection protocol for expression of your target protein
- Uses suspension culture to easily scale up to large amounts of culture
- All reagents are completely animal-origin free, including the defined, serum-free medium, which may be imperative for regulatory requirements.
- Enables transient expression of protein for up to 9 days post-transfection. Peak values vary depending on protein expressed.

For more information, visit [www.lifetechnologies.com](http://www.lifetechnologies.com) or call Technical Support (see page 23).

---

# Methods

## Important guidelines

---

### General cell handling

Follow the general guidelines below to grow and maintain FreeStyle™ 293-F cells.

- **All solutions and equipment that come in contact with the cells must be sterile.** Always use proper sterile technique and work in a laminar flow hood.
  - Before starting experiments, be sure to have cells established (at least 5 passages) and also have some frozen stocks on hand. We recommend using early-passage cells for your experiments (<30 passages). Upon receipt of the cells from Thermo Fisher Scientific, grow and freeze multiple vials of the FreeStyle™ 293-F Cells to ensure that you have an adequate supply of early-passage cells.
  - For general maintenance of cells, pass FreeStyle™ 293-F cells when they reach a density in between  $1 \times 10^6$ – $3 \times 10^6$  viable cells/mL (generally every 48–72 hours). Do not dilute below  $0.1 \times 10^6$  viable cells/mL.
  - Use trypan blue exclusion to determine cell viability (see **Determine cell density and viability**, page 9). Log phase cultures should be >90% viable.
  - When thawing or subculturing cells, transfer cells into pre-warmed medium.
- 

### Important

It is very important to have healthy, well-growing FreeStyle™ 293-F cells to get high yields of protein expression. Strictly follow the guidelines for culturing FreeStyle™ 293-F cells in this manual for the best results.

---



As with other human cell lines, when working with FreeStyle™ 293-F cells, handle as potentially biohazardous material under at least Biosafety Level 2 containment.

---



**Media preparation**

For suspension growth and transfection applications, use:

- FreeStyle™ 293 Expression Medium as is. Supplementation is not required.
  - Antibiotics are **not** recommended; however, 5 mL/L of Antibiotic-Antimycotic containing penicillin, streptomycin, and amphotericin B may be used if required (see page 22 for ordering information).
- 

**Important**

FreeStyle™ 293 Expression Medium is extremely sensitive to light. For optimal results, use and store media protected from light.

---

**Determine cell density and viability**

Follow the procedure below to determine viable and total cell counts.

1. Transfer a small aliquot of the cell suspension to a microcentrifuge tube.
  2. Determine viability and the amount of cell clumping using the trypan blue dye exclusion method (see page 22 for ordering information).
  3. Vigorously vortex for 10–30 seconds to break up cell clumps.
  4. Determine cell density electronically using a Coulter Counter or manually using a hemacytometer.
-

# Thaw and establish cells

---

## Introduction

Follow the protocol below to thaw FreeStyle™ 293-F cells to initiate cell culture. The FreeStyle™ 293-F cell line is supplied in a vial containing 1 mL of cells at  $1 \times 10^7$  viable cells/mL in 90% FreeStyle™ 293 Expression Medium and 10% DMSO. Thaw FreeStyle™ 293-F cells directly into the FreeStyle™ 293 Expression Medium (see page 22 for ordering information).

---

## Materials needed

- FreeStyle™ 293-F cells (supplied with the kit; store frozen cells in liquid nitrogen until ready to use)
  - FreeStyle™ 293 Expression Medium (supplied with the kit; pre-warm to 37°C before use)  
**Note:** We do not recommend adding antibiotics to media as this may negatively impact cell growth.
  - 125-mL polycarbonate, disposable, sterile Erlenmeyer flask with vented cap (available from VWR, Radnor PA, Cat. no. 30180-036)
  - Orbital shaker in 37°C incubator with a humidified atmosphere of 8% CO<sub>2</sub>
  - Reagents to determine viable and total cell counts (see **Determine cell density and viability**, page 9)
-

## Thaw procedure

Store frozen cells in liquid nitrogen until ready to use. To thaw and establish cells:

1. Remove the cryovial of cells from the liquid nitrogen and thaw quickly in a 37°C water bath.
2. Just before the cells are completely thawed, decontaminate the outside of the vial with 70% ethanol. Gently break up clumps and cell pellet if present and transfer the entire contents of the cryovial into a 125-mL polycarbonate, disposable, sterile Erlenmeyer shaker flask containing 30 mL of pre-warmed FreeStyle™ 293 Expression Medium.
3. Incubate cells in a 37°C incubator containing a humidified atmosphere of 8% CO<sub>2</sub> in air on an orbital shaker platform rotating at 135 rpm.
4. Next day, determine viable and total cell counts (see protocol on page 9). Generally, viability is >70%; a bit lower is no reason for concern, but if viability is less than 60% thaw a new batch of cells.
5. Subculture the FreeStyle™ 293-F cells 24–48 hours after thawing by seeding shaker flasks at  $0.3 \times 10^6$  viable cells/mL in pre-warmed FreeStyle™ 293 Expression Medium. We generally use 125- or 250-mL polycarbonate, disposable, sterile, Erlenmeyer flasks containing 40 or 80 mL total working volume of cell suspension, respectively.

**Important Note:** Subculture cells a minimum of 5 passages before use in transfection experiments to allow opportunity for recovery from thawing. To subculture cells, see *Passaging Cells*, page 12.

---

# Subculture cells

---

## Passage cells

Subculture cells when the density is in between  $1-3 \times 10^6$  viable cells/mL, typically every 48–72 hours. When maintaining FreeStyle™ 293-F cells, we generally use a 125- or 250-mL polycarbonate, disposable, sterile Erlenmeyer flask with vented cap containing 25–40 mL or 50–80 mL total working volume of cell suspension, respectively.

**Note:** Glass flasks without baffles may be used, but thorough cleaning after each use is essential to avoid potential toxicity which is more problematic in serum-free cultures.

1. Determine viable and total cell counts (see protocol on page 9).
2. Using the cell density determined in Step 1, calculate the split ratio needed to seed the new shaker flask at  $0.1 \times 10^6$ – $0.2 \times 10^6$  viable cells/mL.
3. Dilute the cells in fresh, pre-warmed FreeStyle™ 293 Expression Medium to give a final cell density of  $0.1 \times 10^6$ – $0.2 \times 10^6$  viable cells/mL in the desired final volume.
4. Incubate flasks in a 37°C incubator containing a humidified atmosphere of 8% CO<sub>2</sub> in air on an orbital shaker platform rotating at 135 rpm.
5. Repeat Steps 1–4 as necessary to maintain or expand cells.



---

FreeStyle™ 293-F suspension cultures may grow as 2–10 cell clusters. Vigorous vortexing for 10–30 seconds may be required at each subculture for a number of passages until the cultures grow predominantly as single cells.

---

## Shake flasks

The cells can be grown in many different culture volumes. We generally use the following polycarbonate, disposable, sterile Erlenmeyer flask with vented cap (other flasks with the same characteristics may be used). For culture volumes **above 400 mL**, lower the speed of the orbital shaker if foam is generated. In 1 L cultures, we recommend 90 rpm.

| Flask Volume | Culture Volume | Manufacturer       | Catalog No. |
|--------------|----------------|--------------------|-------------|
| 125-mL       | 25–40 mL       | VWR, Radnor, PA    | 30180-036   |
| 250-mL       | 50–80 mL       | VWR, Radnor, PA    | 30180-044   |
| 500-mL       | 100–200 mL     | VWR, Radnor, PA    | 30180-052   |
| 1-L          | 200–400 mL     | VWR, Radnor, PA    | 82013-164   |
| 3-L          | 600–1000 mL    | Corning, Acton, MA | 431252      |

**Note:** Glass flasks without baffles may be used, but thorough cleaning after each use is essential to avoid potential toxicity which is more problematic in serum-free cultures.

## Other cell culture systems

It is possible to scale up the FreeStyle™ 293-F cultures in spinner flasks or bioreactors. The appropriate spinner or impeller speed and seeding density should be determined and optimized for each system. At Thermo Fisher Scientific, the optimum spinner speed was 100–130 rpm and 70–100 rpm impeller speed in Celligen™ stirred tank bioreactors. We recommend seeding cells at  $0.3 \times 10^6$ – $0.5 \times 10^6$  viable cells/mL.

**Note:** If the split ratio of cells to fresh media is less than 1:2, you may want to spin down the cell suspension and resuspend the cell pellet in fresh, pre-warmed FreeStyle™ 293 Expression Medium prior to inoculating the spinner or bioreactor culture. Monitor cell viability and the degree of cell clumping. Note that extensive cell clumping may reduce transfection efficiency.



At high stirring speeds (i.e., greater than 130 rpm) and/or depending on the impeller design, you may want to supplement the FreeStyle™ 293 Expression Medium with additional Pluronic™ F-68 (2.5–5 mL/L of 10% Pluronic™ F-68, Cat. no. 24040) to avoid shear stress in the culture.

# Cryopreservation

---

## Introduction

You may freeze FreeStyle™ 293-F cells directly in FreeStyle™ 293 Expression Medium (50:50 ratio of fresh to conditioned medium) with 10% DMSO. When freezing the FreeStyle™ 293-F cell line, we recommend the following:

- Freeze cells at a density of  $1 \times 10^7$  viable cells/mL.
- Use a freezing medium composed of 45% fresh FreeStyle™ 293 Expression Medium, 45% conditioned FreeStyle™ 293 Expression Medium and 10% DMSO.

Guidelines to prepare freezing medium and to freeze cells are provided in this section.

---

## Prepare freezing medium

Prepare freezing medium immediately before use.

1. In a sterile, conical centrifuge tube, mix together the following reagents for every 1 mL of freezing medium needed:

FreeStyle™ 293 Expression Medium

Fresh                      0.45 mL

Conditioned              0.45 mL

DMSO                      0.1 mL

2. Filter-sterilize the freezing medium and place the tube on ice until use. Discard any remaining freezing medium after use.
-

## Freeze cells

Before starting, label cryovials and prepare freezing medium. Keep the freezing medium on ice.

1. Grow the desired quantity of FreeStyle™ 293-F cells in shaker flasks, harvesting when the cell density reaches  $0.5 \times 10^6$ – $1 \times 10^6$  viable cells/mL. Transfer cells to a sterile, conical centrifuge tube.
2. Determine the viable and total cell counts (see protocol on page 9) and calculate the volume of freezing medium required to yield a final cell density of  $1 \times 10^7$  viable cells/mL.
3. Centrifuge cells at  $100 \times g$  for 5 minutes at room temperature and carefully aspirate the medium.
4. Resuspend the cells in the pre-determined volume of chilled freezing medium.
5. Place cryovials in a microcentrifuge rack and aliquot 1 mL of the cell suspension into each cryovial.
6. Freeze cells in an automated or manual, controlled-rate freezing apparatus following standard procedures. For ideal cryopreservation, the freezing rate should be a decrease of  $1^\circ\text{C}$  per minute.
7. Transfer frozen vials to liquid nitrogen for long-term storage.

**Note:** You may check the viability and recovery of frozen cells 24 hours after storing cryovials in liquid nitrogen by following the procedure outlined in **Thaw and establish cells**, page 10.

---

# Transfect cells

---

## Introduction

To transfect suspension FreeStyle™ 293-F cells, you will use the cationic lipid-based transfection reagent, FreeStyle™ MAX Reagent, complex in animal-origin free OptiPro™ SFM, while cells stay in FreeStyle™ 293 Expression Medium, which does not inhibit cationic lipid-mediated transfection. FreeStyle™ 293 Expression Medium is formulated to allow high transfection efficiency of suspension FreeStyle™ 293-F cells **without** the need to change media. Transient transfection experiments may be performed in a large volume, allowing large-scale protein production.

FreeStyle™ 293 Expression Medium, FreeStyle™ MAX Reagent and OptiPro™ SFM are available separately from Thermo Fisher Scientific (see page 22 for ordering information). For more information, see the FreeStyle™ MAX 293 Expression System, available at [www.lifetechnologies.com](http://www.lifetechnologies.com) or call Technical Support (see page 23).

---

## FreeStyle™ MAX Reagent

FreeStyle™ MAX Reagent is a proprietary formulation suitable for transfection of DNA into eukaryotic cells grown in suspension and provides the following advantages:

- FreeStyle™ MAX Reagent demonstrates high transfection efficiency with minimal cytotoxicity in suspension FreeStyle™ 293-F cells (cultured in FreeStyle™ 293 Expression Medium)
  - DNA-FreeStyle™ MAX Reagent complexes can be added directly to cells in culture medium
  - It is not necessary to remove complexes or change or add medium following transfection
- 

## Plasmid preparation

Plasmid DNA for transfection into eukaryotic cells must be clean, sterile and free from contamination with phenol and NaCl. Contaminants may kill the cells, and salt will interfere with complexing, decreasing transfection efficiency. We recommend isolating DNA using the PureLink™ HiPure Plasmid Kits, which are validated for use with the FreeStyle™ MAX 293 Expression System (see page 22).

**Note:** Make sure your DNA preparation is sterile. We recommend performing filtration before use through a 0.22-µm filter.

---



## Materials needed

- Suspension FreeStyle™ 293-F cells cultured in FreeStyle™ 293 Expression Medium.  
**Recommendation:** Calculate the number of cells that you will need for your transfection experiment and expand cells accordingly. Make sure that the cells are healthy and greater than 95% viable before proceeding to transfection.
- Purified plasmid DNA of interest (1 mg/mL).
- FreeStyle™ MAX Reagent (store at 4°C until use).
- OptiPro™ SFM to dilute DNA and transfection reagent (pre-warmed to room temperature). See page 22 for ordering information.
- FreeStyle™ 293 Expression Medium (pre-warmed to 37°C).  
**Note:** Do not add antibiotics to media during transfection as this may decrease transfection activity.
- 125-mL polycarbonate, disposable, sterile Erlenmeyer flasks.
- Orbital shaker in 37°C incubator with a humidified atmosphere of 8% CO<sub>2</sub>.
- Reagents to determine viable and total cell counts.

FreeStyle™ 293 Expression Medium, OptiPro™ SFM and FreeStyle™ MAX Reagent are available separately from Thermo Fisher Scientific (see page 22 for ordering information).

---

## Optimal conditions for 30 mL transfection

To transfect suspension FreeStyle™ 293-F cells in a 30 mL volume, we recommend using the following optimized conditions:

- **Final transfection volume:** 30 mL
  - **Number of cells to transfect:**  $3 \times 10^7$  cells (final cell density of  $1 \times 10^6$  cells/mL)
  - **Amount of plasmid DNA:** 37.5 µg (starting point; can vary from 24–42 µg)
  - **FreeStyle™ MAX Reagent:** 37.5 µL (starting point; can vary from 24–42 µL)
-

## Transfection procedure

Follow the procedure below to transfect suspension FreeStyle™ 293-F cells in a 30 mL volume. Remember that you may keep the cells in FreeStyle™ 293 Expression Medium during transfection. We recommend including a positive control (pCMV SPORT-βgal) and a negative control (no DNA, no FreeStyle™ MAX Reagent) in your experiment to help you evaluate your results.

1. Approximately 24 hrs before transfection, pass FreeStyle™ 293-F cells at  $0.6 \times 10^6$ – $0.7 \times 10^6$  cells/mL. Place the flask(s) on an orbital shaker platform rotating at 135 rpm at 37°C, 8% CO<sub>2</sub>.
  2. On the day of transfection, the cell density should be about  $1.2 \times 10^6$ – $1.5 \times 10^6$ /mL. Dilute the cells to  $1 \times 10^6$  cells/mL. To ensure high transfection results, viability of cells must be over 90%. Add 30 mL of cells into each 125-mL shake flask.
  3. Gently invert the tube of FreeStyle™ MAX Transfection Reagent several times to mix. Do not vortex.
  4. Dilute 37.5 μg of plasmid DNA into OptiPro™ SFM to a total volume of 0.6 mL and mix. In a separate tube, dilute 37.5 μL of FreeStyle™ MAX Reagent in OptiPro™ SFM to a total volume of 0.6 mL and mix gently by inverting the tube (do not vortex). Immediately add diluted FreeStyle™ MAX Reagent to diluted DNA solution to obtain a total volume of 1.2 mL and mix gently.
  5. Incubate the DNA-lipid mixture for 10 minutes at room temperature to allow complexes to form. Do not incubate for longer than 20 minutes.
  6. Slowly add 1.2 mL of DNA-lipid mixture into the 125-mL flask containing cells while slowly swirling the flask.
  7. Incubate transfected cell cultures at 37°C, 8% CO<sub>2</sub> on an orbital shaker platform rotating at 135 rpm. There is no need to change or supplement the culture medium during the first 6–7 days.
  8. Protein expression may be detectable within 4–8 hours of transfection, with maximal protein yield usually between 1–7 days post-transfection, depending on the protein expressed.
-

## Optimize protein expression

- When expressing a protein for the first time, perform a time course experiment between days 1 and 9 post-transfection to identify the peak of protein production, and to monitor cell viability.
- Test varying amounts of plasmid DNA and FreeStyle™ MAX Reagent. For 30 mL cultures, try between 24–42 µg DNA and 24–42 µL FreeStyle™ MAX Reagent.
- For secreted IgG protein production, we have observed peak yields at 5–7 days post-transfection.
- To assess transfection efficiency via expression of a GFP-type fluorescent protein, we recommend monitoring the cultures starting at 24 hours post-transfection.
- In some cases, transfection efficiency may go down during the course of the experiment, although protein production is still going up. Never evaluate results by transfection efficiency only; always measure the protein production.

For optimizing protein expression while scaling up culture volumes, see **Scale up transfections** in the following section.

**Note:** Cells that are transfected with high efficiency grow slower than untransfected cells. This is a good thing.

---

## Scale up transfections

If you transfect suspension FreeStyle™ 293-F cells in a larger volume, consider the following points:

- Scale up the volume of each reagent in proportion to the culture volume.
  - For culture volumes **above 40 mL**, lower the speed of the orbital shaker if foam is generated. In 1 L cultures, we recommend 90 rpm.
  - The transfection conditions may vary depending on the type of culture vessel used and the growth conditions of your cells; therefore, you may want to perform pilot studies to optimize your transfection conditions.
  - Often, the conditions for transfection of large culture volumes need to be adjusted to the higher side of the range, while those for small culture volumes may tend towards the lower side of the range.
-

# Troubleshooting

**Culture cells** The following table lists some potential problems and solutions that help you troubleshoot your cell culture problems.

| Problem  | Cause   | Solution  |
|--|---|---|
| No viable cells after thawing stock  | Stock not stored correctly                                    | Order new stock and store in liquid nitrogen. Keep in liquid nitrogen until thawing.  |
|  | Home-made stock not viable                                    | Freeze cells at a density of $1 \times 10^7$ viable cells/mL.   |
|  |   | Use a freezing medium composed of 45% fresh FreeStyle™ 293 Expression Medium, 45% conditioned FreeStyle™ 293 Expression Medium and 10% DMSO.          |
|  |   | Use low-passage cells to make your own stocks.  |
|  |   | Follow procedures in <b>Cryopreservation</b> (page 14).   |
|  | Obtain new FreeStyle™ 293-F Cells.                            |   |
|  | Thawing medium not correct                                    | Use FreeStyle™ 293 Expression Medium (pre-warm before use)  |
| Do not add antibiotics to media as this may negatively impact cell growth. |   |   |
| Shaker not set up properly   | See page 12–13 for proper settings for orbital shaker.        |   |
| Cells too diluted  | Spin down culture and grow cells in a smaller culture volume. |   |
| Cells grow slowly  | Growth medium not correct                                     | Use FreeStyle™ 293 Expression Medium (pre-warm before use).   |
|  |   | Do not add antibiotics to media as this may negatively impact cell growth.  |
|  | Shaker not set up properly                                    | See page 12 for proper settings for orbital shaker.   |
|  | Medium foamy  | Lower the shaker speed slightly till no foam forms.   |
|  | Flasks too small  | Use flasks that are at least 2.5 times bigger than the culture volume.  |
|  | Cells too old   | Use healthy FreeStyle™ 293-F cells under passage 30; do not overgrow.   |
|  | Cell culture clumpy   | Prevent this by sufficient agitation of the culture, a regular and frequent cell passage schedule, and maintenance of cells at recommended densities. |
|  | Cell density too low when passing cells                       | Spin down cells at $100 \times g$ for 5 minutes at room temperature, aspirate media and dilute cells at higher density.                               |
|  | Cells transfected   | Cells transfected at high efficiency grow slower than untransfected cells. This is a desired outcome.   |

## Transfection and protein production

The following table lists some potential problems and possible solutions that may help you troubleshoot your transfection and protein production experiments.

| Problem  | Cause  | Solution   |  |
|--|--|--|--|
| Low Transfection Efficiency and/or Low Protein Yield | Cells cultured for too many passages ( $\geq 30$ ) | Thaw a new batch of early-passage cells.   |  |
|  | Cells not passed 24 hours before transfection      | Approximately 24 hours before transfection, pass cells at $6\text{--}7 \times 10^5$ cells/mL.  |  |
|  | Improperly cultured FreeStyle™ 293-F cells         | Exactly follow procedures as outlined in <b>Subculture Cells</b> section (page 12).  |  |
|  | Cells transfected in media containing antibiotics  | Do not add antibiotics during transfection.  |  |
|  | FreeStyle™ Max Reagent handled incorrectly         |  | Store at 4°C. <b>Do not freeze.</b>  |
|  |  |  | Mix gently by inversion. <b>Do not vortex.</b>   |
|  | Used poor quality expression construct plasmid DNA | <b>Do not</b> use mini-prep plasmid DNA for transfection. Use a PureLink™ HiPure Plasmid Kit to prepare plasmid DNA with low endotoxin contamination.  |  |
|  | Suboptimal transfection conditions                 |  | Perform transfections with positive control plasmid pCMV SPORTβ-□gal to <b>assess</b> your transfection conditions.  |
|  |  |  | Assess transfection efficiency via expression of a GFP-type fluorescent protein (we recommend monitoring the cultures starting at 24 hours post-transfection). |
|  |  |  | Vary the amounts of DNA and FreeStyle™ MAX Reagent used (see page 19)  |
| DNA not sterile                                      |  | Sterilize DNA (see page 16).   |  |
| Gene of interest is toxic to cells                   |  | Do not generate constructs containing activated oncogenes or harmful genes.  |  |
|  |  | Try FreeStyle™ MAX CHO Expression System.  |  |
| Protein harvested too early or too late              |  | When expressing a protein for the first time, perform a time course experiment between days 1 and 7 post-transfection to identify the peak of protein production, and to monitor cell viability. |  |

# Appendix

## Accessory products

### Additional products

The products listed in this section may be used with the FreeStyle™ 293-F Cells. For more information, refer to our website ([www.lifetechnologies.com](http://www.lifetechnologies.com)) or call Technical Support (see page 23).

| Item  | Quantity    | Catalog no. |
|---|-------------|-------------|
| FreeStyle™ MAX 293 Expression System                      | 1 kit       | K9000-10    |
| FreeStyle™ 293 Expression Medium                          | 1 L         | 12338-018   |
|   | 6 × 1 L     | 12338-026   |
| FreeStyle™ MAX Reagent                                    | 1 mL        | 16447-100   |
| OptiPro™ SFM  | 100 mL      | 12309-050   |
|   | 1000 mL     | 12309-019   |
| PureLink™ HiPure Plasmid Midiprep Kit                     | 25 preps    | K2100-04    |
| PureLink™ HiPure Plasmid Maxiprep Kit                     | 10 preps    | K2100-06    |
| PureLink™ HiPure Plasmid Filter Maxiprep Kit              | 10 preps    | K2100-16    |
| PureLink™ HiPure Plasmid Megaprep Kit                     | 4 preps     | K2100-08    |
| Trypan Blue Stain   | 100 mL      | 15250-061   |
| FluoReporter™ <i>lacZ</i> /Galactosidase Quantitation Kit | 1000 assays | F-2905      |
| Antibiotic-Antimycotic (100X)                             | 100 mL      | 15240-062   |

### FreeStyle™ MAX CHO Expression System

Protein production using the FreeStyle™ MAX Expression Systems can be performed in CHO cells or 293 cells. The following table lists ordering information for the reagents specific to the FreeStyle™ MAX CHO Expression System. Note that the FreeStyle™ MAX Reagent is used with both the FreeStyle MAX 293 and CHO Systems.

| Item                                 | Quantity                           | Catalog no. |
|--------------------------------------|------------------------------------|-------------|
| FreeStyle™ MAX CHO Expression System | 1 kit                              | K9000-20    |
| FreeStyle™ CHO-S Cells               | 1 vial (1 × 10 <sup>7</sup> cells) | R800-07     |
| FreeStyle™ CHO Expression Medium     | 1 L                                | 12651-014   |
|                                      | 6 × 1 L                            | 12651-022   |

# Technical support

---

## Obtaining support

For the latest services and support information for all locations, go to **[www.lifetechnologies.com](http://www.lifetechnologies.com)**.

At the website, you can:

- Access worldwide telephone and fax numbers to contact Technical Support and Sales facilities
  - Search through frequently asked questions (FAQs)
  - Submit a question directly to Technical Support (**[techsupport@lifetech.com](mailto:techsupport@lifetech.com)**)
  - Search for user documents, SDSs, vector maps and sequences, application notes, formulations, handbooks, certificates of analysis, citations, and other product support documents
  - Obtain information about customer training
  - Download software updates and patches
- 

## Safety Data Sheets (SDS)

Safety Data Sheets (SDSs) are available at **[www.lifetechnologies.com/support](http://www.lifetechnologies.com/support)**.

---

## Certificate of analysis

The Certificate of Analysis provides detailed quality control and product qualification information for each product. Certificates of Analysis are available on our website. Go to **[www.lifetechnologies.com/support](http://www.lifetechnologies.com/support)** and search for the Certificate of Analysis by product lot number, which is printed on the box.

---

## Limited product warranty

Life Technologies and/or its affiliate(s) warrant their products as set forth in the Life Technologies General Terms and Conditions of Sale found on the Life Technologies web site at **[www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions)**. If you have any questions, please contact Life Technologies at **[www.lifetechnologies.com/support](http://www.lifetechnologies.com/support)**.

---

# Purchaser notification

---

## Information for European customers

The FreeStyle™ 293-F Cell Line is genetically modified and carries human adenovirus type 5 DNA. As a condition of sale, this product must be in accordance with all applicable local legislation and guidelines including EC Directive 90/219/EEC on the contained use of genetically modified organisms.

---



# References

---

- Graham, F. L., Smiley, J., Russell, W. C., and Nairn, R. (1977) Characteristics of a Human Cell Line Transformed by DNA from Human Adenovirus Type 5. *J. Gen. Virol.* 36, 59-74
- Harrison, T., Graham, F., and Williams, J. (1977) Host-range Mutants of Adenovirus Type 5 Defective for Growth in HeLa Cells. *Virology* 77, 319-329
-

For support visit [lifetechnologies.com/support](https://lifetechnologies.com/support)  
or email [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

[lifetechnologies.com](https://lifetechnologies.com)

**ThermoFisher**  
SCIENTIFIC