


# Poly(A) Polymerase (Cloned)

Catalog Number AM2030

Pub. No. 4393885 Rev. C

Contents	Quantity	Storage conditions
Poly(A) Polymerase, 2 U/μL	80 Units	Store at -20°C. Do not store in a frost-free freezer.
5X E-PAP Buffer	800 μL	
25 mM MnCl <sub>2</sub>	400 μL	
DEPC treated Water	1 mL	

 **WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## Product description

Poly(A) Polymerase (PAP) catalyzes the addition of adenosine to the 3' end of RNA in a sequence-independent fashion.

**Source:** An *E. coli* strain overexpressing *E. coli* Poly(A) Polymerase I.

**Unit (U) definition:** One unit is the amount of enzyme that incorporates 1 nmol of AMP into an acid insoluble tail on tRNA in 10 min at 37°C at pH 7.9 with ATP as the substrate.

**Storage buffer (not included):** 25 mM Tris (pH 8.0), 500 mM NaCl, 0.1 mM DTT, 1 mM MgCl<sub>2</sub>, 0.1 mM EDTA and 50% glycerol (v/v).

## Using Poly(A) Polymerase

Poly(A) Polymerase can be used to add a poly(A) tail to RNA transcripts. A poly(A) tail may increase mRNA stability and therefore translation efficiency when tailed RNA is used in transfection or micro-injection experiments. It can also be used to label the 3' end of RNA molecules by addition of radiolabeled adenosine.

### Poly(A) tailing reactions

Use 8 U Poly(A) Polymerase to generate ≥150 nt-long poly(A) tails. Use less Poly(A) Polymerase to produce shorter poly(A) tails.

In a total reaction volume of 100 μL:

- ≤40 μg RNA
- 1X PAP Buffer
- 2.5 mM MnCl<sub>2</sub>
- 1 mM ATP
- 2–8 U Poly(A) Polymerase

For additional information on poly(A) tailing reactions, refer to the Poly(A) Tailing Kit (Cat. no. AM1350) user guide, available at [www.lifetechnologies.com](http://www.lifetechnologies.com).

### Other applications

PAP can be used to append radiolabeled adenosine to the 3' end of RNA molecules to generate labeled RNAs.

In conjunction with terminal transferase, reverse transcriptase, and thermostable polymerase, PAP can be used to clone RNAs of unknown sequence (see the following figure). For RNA cloning, first add a poly(A) tail to the 3' end of the RNA using PAP. Reverse transcribe the RNA using reverse transcriptase (for example, Cat. no. AM2043 or AM2044) and an oligo(dT) primer. Add a poly(dA) tail to the 3' end of the cDNA using terminal transferase. Then amplify the modified cDNA using SuperTaq™ DNA Polymerase (Cat. no. AM2052, AM2054)

and PCR primers specific to the homopolymer tails appended by PAP and terminal transferase. The dsDNA can then be cloned and sequenced. Using PCR primers with restriction sites can facilitate cloning.

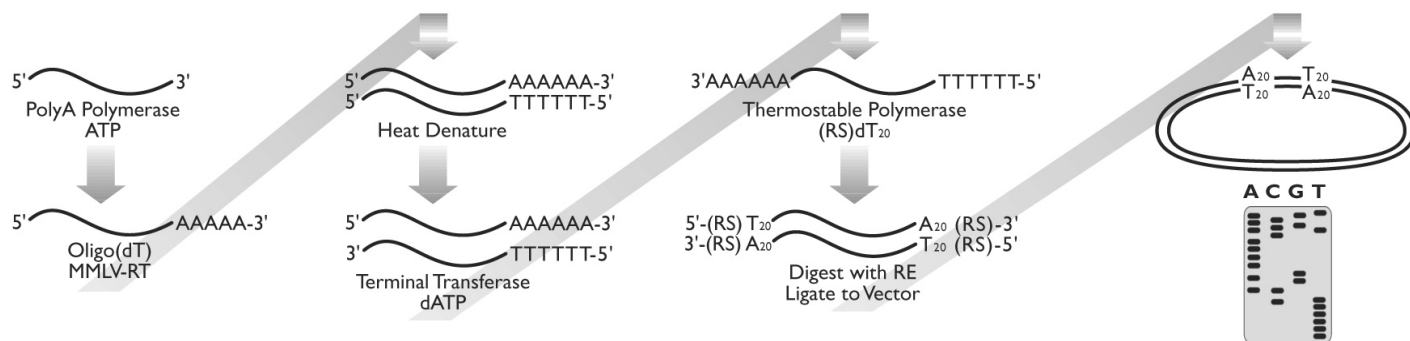


Figure 1 Cloning RNAs with PAP

## Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at [www.lifetechnologies.com/termsandconditions](http://www.lifetechnologies.com/termsandconditions). If you have any questions, please contact Life Technologies at [www.lifetechnologies.com/support](http://www.lifetechnologies.com/support).

## References

Biebricher, C.K. and Luce, R. 1993. Sequence analysis of RNA species synthesized by Q beta replicase without template. *Biochemistry* 32: 4848–4854.

### For Research Use Only. Not for use in diagnostic procedures.

The information in this guide is subject to change without notice.

DISCLAIMER: LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

NOTICE TO PURCHASER: LIMITED USE LABEL LICENSE: Research Use Only : The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact [outlicensing@lifetech.com](mailto:outlicensing@lifetech.com) or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

© 2012 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners. SuperTaq is a trademark of Enzyme Technologies Ltd.

For support visit [lifetechnologies.com/support](http://lifetechnologies.com/support) or email [techsupport@lifetech.com](mailto:techsupport@lifetech.com)

lifetechnologies.com

2 November 2012

