CellTrace™ Violet Cell Proliferation Kit

Cat. no. C34557, C34571

Pub no MANDON6888

Revision A.0

Detailed protocol is available online at www.thermofisher.com.

WARNING! For every chemical, read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from www.thermofisher.com/support.

Chemical	CAS	Warning
DMS0	67-68-5	Components of the product may be absorbed into the body through the skin.



1. Remove supernatant from the cell pellet.



2. Add CellTrace™ Violet (1:1000 dilution) staining solution and gently re-suspend the cells.



3. Incubate at 37°C for 20 minutes, protected from light.



4. Add complete culture medium and mix.



5. Incubate at 37°C for 5 minutes.



6. Pellet the cells and remove the supernatant.



7. Resuspend the cell pellet in fresh, pre-warmed complete culture medium.



📕 🔄 8. Analyze by flow cytometry or proceed with stimulation or incubation

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support

DISCLAIMER - TO THE EXTENT ALLOWED BY LAW LIFE TECHNOLOGIES AND/OR ITS AFEILIATE(S) WILL NOT BE LIABLE FOR SPECIAL INCIDENTAL INDIRECT. PLINITIVE MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

© 2015 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

For support visit thermofisher.com/support