## ProLong® Gold Antifade Reagent ProLong® Gold Antifade Reagent with DAPI

Pub. no. MAN0009669

Revision A

Detailed protocol is available online at www.lifetechnologies.com/manuals.

4 0 1		
1 × 2 mL	P10144	Room temperature (15–30°C)     Protect from light
5 × 2 mL	P36934	
× 10 mL	P36930	
1 × 2 mL	P36941	
5 × 2 mL	P36935	
× 10 mL	P36931	
5	× 10 mL × 2 mL 5 × 2 mL	× 10 mL P36930 × 2 mL P36941 5 × 2 mL P36935 × 10 mL P36931

<sup>\*</sup> Product may also be stored at <-20°C.

life technologies"

For Research Use Only. Not for use in diagnostic procedures.



1. Add 1 drop of ProLong® Gold reagent onto glass slide.



2. Remove excess liquid from coverslip.



3. Place coverslip sample-side down onto ProLong® Gold reagent on the glass slide.



4. Cure for 24 hours at room temperature in the dark



5. Clean coverslip and image cells.

## Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support. DISCLAIMER - LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATEISI DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT. EXPRESSED OR IMPLIED. INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTARILITY FITNESS FOR A PARTICULAR PURPOSE OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATEISLE LIARLE WHETHER IN CONTRACT TORT WARRANTY OR LINDER ANY STATLITE OR ON ANY OTHER BASIS FOR SPECIAL INCIDENTAL INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF © 2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or

For support visit lifetechnologies.com/support