



# **Labeled Goat Anti-Mouse IgM Antibodies**

**Table 1.** Contents and storage information.

Material	Amount	Concentration	Storage	Stability *
Alexa Fluor® dye and DSB-X™ biotin conjugates	250 μL	2 mg/mL in 0.1 M sodium phosphate, 0.1 M NaCl, pH 7.5, 5 mM sodium azide	Store undiluted 2–6°C Protect from light Avoid repeated freezing and thawing	When stored as directed, products are stable for at least 3 months.
Allophycocyanin (APC) and R-phycoerythrin (RPE) conjugates		1 mg/mL in 0.1 M sodium phosphate, 0.1 M NaCl, pH 7.5, 5 mM sodium azide		

<sup>\*</sup> For longer storage, divide the solution into single-use aliquots and freeze at ≤-20°C. Frozen aliquots are stable for at least six months. AVOID REPEATED FREEZING AND THAWING.

Degree of labeling: Typically 3–8 fluorophore or biotin molecules per antibody; the exact degree of labeling is indicated on the product

Approximate fluorescence excitation/emission maxima: See Table 2.

## Introduction

Invitrogen's Molecular Probes™ goat anti-mouse IgM conjugates (Table 2) are prepared from antibodies purified by mouse IgM-affinity chromatography. The antibodies react specifically with IgM heavy chains (µ chains) and not with immunoglobulin light chains. To minimize cross-reactivity, the goat anti-mouse IgM antibodies have been adsorbed against human IgM prior to labeling.

In addition to the antibodies listed in this product manual, Invitrogen prepares conjugates of many species-specific anti-IgG antibodies, as well as conjugates of avidin, streptavidin, NeutrAvidin™ biotin-binding protein, protein A, and protein G. Refer to www.invitrogen.com or contact Technical Support for more information about these products.

At the time of preparation, the products are certified to be free of unconjugated label.

Table 2. Invitrogen's labeled goat anti-mouse IgM antibodies.\*

Catalog #	Label	Ex †	Em †
A31552	Alexa Fluor® 350	346	442
A21042	Alexa Fluor® 488	495	519
A21045	Alexa Fluor® 546	556	573
A21426	Alexa Fluor® 555	555	565
A21043	Alexa Fluor® 568	578	603
A21044	Alexa Fluor® 594	590	617
A21046	Alexa Fluor® 633 ‡	632	647
A21238	Alexa Fluor® 647 ‡	650	668
A21048	Alexa Fluor® 680 ‡	679	702
A10688	allophycocyanin	650	660
A10689	R-phycoerythrin (RPE)	496	578
A10690 §	R-phycoerythrin (RPE)	496	578
D20693	DSB-X™ biotin	NA	NA

<sup>\*</sup>These goat anti-mouse IgM antibodies have been adsorbed against human  $IgG_1$ ,  $IgG_{2a}$ ,  $IgG_{2b}$ ,  $IgG_3$ , IgA, human serum and purified human para-proteins to minimize cross-reactivity. † Approximate fluorescence excitation (Ex) and emission (Em) maxima, in nm. ‡ Human vision is insensitive to light beyond ~650 nm, and therefore it is not possible to view the fluorescence of the Alexa Fluor® 633, 647, and 680 dyes by looking through a conventional fluorescence microscope. § F(ab')2 fragment.

## **Guidelines for Use**

It is a good practice to centrifuge the protein conjugate solution briefly in a microcentrifuge before use; only the supernatant should then be added to the experiment. This step eliminates any protein aggregates that may form during storage, and reduces nonspecific background staining.

Because staining protocols vary with application, the appropriate dilution of antibody should be determined empirically. For these labeled antibodies, a final concentration of  $1-10~\mu g/mL$ should be satisfactory for most immunohistochemical applications.<sup>1</sup>

## Reference

1. Short Protocols in Molecular Biology, 2nd Edition, F.M. Ausubel et al., Eds., John Wiley and Sons (1992) pp. 14-24–14-30.

## Product List Current prices may be obtained from our website or from our Customer Service Department.

Cat #	Product Name	Unit Size
A31552	Alexa Fluor® 350 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21042	Alexa Fluor® 488 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21045	Alexa Fluor® 546 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21426	Alexa Fluor® 555 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21043	Alexa Fluor® 568 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21044	Alexa Fluor® 594 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21046	Alexa Fluor® 633 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21238	Alexa Fluor® 647 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
A21048	Alexa Fluor® 680 goat anti-mouse IgM (μ chain) *2 mg/mL*	250 μL
D20693	DSB-X™ biotin goat anti-mouse lgM (µ chain) *2 mg/mL*	250 μL

## **Contact Information**

#### Molecular Probes, Inc.

29851 Willow Creek Road Eugene, OR 97402 Phone: (541) 465-8300 Fax: (541) 335-0504

#### **Customer Service:**

6:00 am to 4:30 pm (Pacific Time) Phone: (541) 335-0338 Fax: (541) 335-0305 probesorder@invitrogen.com

#### **Toll-Free Ordering for USA:**

Order Phone: (800) 438-2209 Order Fax: (800) 438-0228

#### **Technical Service:**

8:00 am to 4:00 pm (Pacific Time) Phone: (541) 335-0353 Toll-Free (800) 438-2209 Fax: (541) 335-0238 probestech@invitrogen.com

#### **Invitrogen European Headquarters**

Invitrogen, Ltd. 3 Fountain Drive Inchinnan Business Park Paisley PA4 9RF, UK Phone: +44 (0) 141 814 6100 Fax: +44 (0) 141 814 6260 Email: euroinfo@invitrogen.com Technical Services: eurotech@invitrogen.com Further information on Molecular Probes products, including product bibliographies, is available from your local distributor or directly from Molecular Probes. Customers in Europe, Africa and the Middle East should contact our office in Paisley, United Kingdom. All others should contact our Technical Service Department in Eugene, Oregon.

Molecular Probes products are high-quality reagents and materials intended for research purposes only. These products must be used by, or directly under the supervision of, a technically qualified individual experienced in handling potentially hazardous chemicals. Please read the Material Safety Data Sheet provided for each product; other regulatory considerations may apply.

### Limited Use Label License No. 223: Labeling and Detection Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (b) its components to a third party or otherwise transfer (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (b) its components or (c) materials made using this product or its components to a third party or otherwise transfer (b) its components or (c) materials made using the components of the component of the componen erwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) to not transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for the rapeutic. diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is a compared to the manufacture of such products of the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is a compared to the manufacture of such products of the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is a compared to the purchaser is not will be used to the purchaser in the purchaser is not will be used to the purchaser in the purchaser is not will be used to the purchaser in the purchaser in the purchaser is not will be used to the purchaser in the purchaser in the purchaser is not will be used to the purchaser in the purchasewilling to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Molecular Probes, Inc., Business Development, 29851 Willow Creek Road, Eugene, OR 97402, Tel: (541) 465-8300. Fax: (541) 335-0354.

Several Molecular Probes products and product applications are covered by U.S. and foreign patents and patents pending. All names containing the designation  $\ensuremath{^\circ}$  are registered with the U.S. Patent and Trademark Office.

Copyright 2009, Molecular Probes, Inc. All rights reserved. This information is subject to change without notice.