



ELISA Kit  
Catalog #KHO0761

**Human  
STAT5a  
[pY694]\***

[www.invitrogen.com](http://www.invitrogen.com)

Invitrogen Corporation

542 Flynn Road, Camarillo, CA 93012

Tel: 800-955-6288

**E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)**

\*Patent Pending



## TABLE OF CONTENTS

Introduction .....	4
Principle of the Method.....	5
Reagents Provided.....	6
Supplies Required But Not Provided .....	7
Procedural Notes/Lab Quality Control.....	7
Safety .....	9
Directions for Washing.....	9
Procedure for Extraction of Proteins from Cells.....	10
Reagent Preparation and Storage.....	12
Reconstitution and Dilution of Hu STAT5a [pY694] Standard.....	12
Dilution of Hu STAT5a [pY694] Standard.....	13
Storage and Final Dilution of Anti-Rabbit IgG HRP (100X).....	14
Dilution of Wash Buffer .....	15
Assay Method.....	15
Typical Data .....	18
Limitations of the Procedure.....	19
Performance Characteristics .....	20
Sensitivity .....	20
Precision .....	21
Recovery.....	22
Parallelism .....	22
Linearity of Dilution .....	23
Specificity.....	24
References .....	27

## INTRODUCTION

Signal Transducer and Activator of Transcription 5 (STAT5, consists of isoforms a and b) is a ~95 kDa member of the STAT family of proteins, which is comprised of STATs 1, 2, 3, 4, 5a, 5b, and 6. STAT homologs have also been described in *Drosophila* sp. and *Dictyostellium discoideum*. STAT family members are characterized by the presence of a DNA binding domain, an SH3 domain which mediates interaction with polyproline-containing proteins, an SH2 domain which mediates interaction with phosphotyrosine-containing proteins, and a C-terminal transactivation (TAD) domain. STAT proteins control the transcription of specific genes in response to cytokine stimulation.

STAT5a was originally identified in lactating mammary gland and is the primary transcription factor responsible for signaling by prolactin (Prl). Following stimulation by Prl, the principal STAT5 isoform, STAT5a, is expressed in the mammary gland and is rapidly tyrosine phosphorylated by the JAK2 kinase, dimerizes and translocates into the nucleus where it interacts as a tetramer with clustered STAT5a binding sites in both the  $\beta$ -casein promoter and enhancer.

STAT5a was subsequently found to have diverse functions in many cell types, where it is activated to regulate gene transcription in response to multiple hematopoietic cytokines and growth factors. STAT5 is involved in normal cell cycle, survival, and apoptosis, and is implicated in leukaemogenesis and angiogenesis. STAT5 is activated by phosphorylation of tyrosine 694 in STAT5a (tyrosine 699 in STAT5b), a major Src phosphorylation site that is critical for STAT5 dimerization, translocation and binding to DNA.

The Invitrogen STAT5a [pY694] Human ELISA is designed to detect and quantify the level of STAT5a protein phosphorylated at tyrosine residue 694. This assay is intended for the detection of STAT5a [pY694] from lysates of human cells or tissues. Reactivity with other species has not been thoroughly evaluated. Invitrogen also offers a STAT5a (Total) ELISA kit (Cat. # KHO0751), which quantifies STAT5a independently of phosphorylation status and allows normalization of phosphorylated STAT5a to total STAT5a.

**For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.**

## **Read entire protocol before use**

### **PRINCIPLE OF THE METHOD**

The Invitrogen STAT5a [pY694] kit is a solid phase sandwich Enzyme Linked-Immuno-Sorbent Assay (ELISA). A monoclonal antibody specific for STAT5a (regardless of phosphorylation state) has been coated onto the wells of the microtiter strips provided. Samples, including a standard containing STAT5a [pY694], control specimens, and unknowns, are pipetted into these wells. During the first incubation, the STAT5a antigen binds to the immobilized (capture) antibody. After washing, a rabbit antibody specific for STAT5a phosphorylated at tyrosine 694 is added to the wells. During the second incubation, this antibody serves as a detection antibody by binding to the immobilized STAT5a [pY694] protein captured during the first incubation. After removal of excess detection antibody, a horseradish peroxidase-labeled Anti-Rabbit IgG (Anti-Rabbit IgG HRP) is added. This binds to the detection antibody to complete the four-member sandwich. After a third incubation and washing to remove all the excess Anti-Rabbit IgG HRP, a substrate solution is added, which is acted upon by the bound enzyme

to produce color. The intensity of this colored product is directly proportional to the concentration of STAT5a [pY694] present in the original specimen.

## REAGENTS PROVIDED

**Note:** Store all reagents at 2 to 8°C.

<i>Reagent</i>	<i>96 Test Kit</i>
<i>Hu STAT5a [pY694] Standard.</i> Contains 0.1% sodium azide. Refer to vial label for quantity and reconstitution volume.	2 vials
<i>Standard Diluent Buffer.</i> Contains 0.1% sodium azide; red dye*; 25 mL per bottle.	1 bottle
<i>Antibody Coated Wells.</i> 12 x 8 Well Strips.	1 plate
<i>Hu STAT5a [pY694] Detection Antibody.</i> Contains 0.1% sodium azide; blue dye*; 11 mL per bottle.	1 bottle
<i>Anti-Rabbit IgG HRP (100X).</i> Contains 3.3 mM thymol; 0.125 mL per vial.	1 vial
<i>HRP Diluent.</i> Contains 3.3 mM thymol; yellow dye*; 25 mL per bottle.	1 bottle
<i>Wash Buffer Concentrate (25X).</i> 100 mL per bottle.	1 bottle
<i>Stabilized Chromogen, Tetramethylbenzidine (TMB).</i> 25 mL per bottle.	1 bottle
<i>Stop Solution.</i> 25 mL per bottle.	1 bottle
<i>Plate Covers,</i> adhesive strips.	3
* In order to help our customers avoid any mistakes in pipetting the ELISAs, we provide colored <i>Standard Diluent Buffer</i> , <i>Detection Antibody</i> , and <i>HRP Diluent</i> to help monitor the addition of solutions to the reaction wells. This does not in any way interfere with the test results.	

**Disposal Note:** This kit contains materials with small quantities of sodium azide. Sodium azide reacts with lead and copper plumbing to form explosive metal azides. Upon disposal, flush drains with a large volume of water to prevent azide accumulation. Avoid ingestion and contact with eyes, skin and mucous membranes. In case of contact, rinse affected area with plenty of water. Observe all federal, state and local regulations for disposal.

### **SUPPLIES REQUIRED BUT NOT PROVIDED**

1. Microtiter plate reader capable of measurement at or near 450 nm.
2. Calibrated adjustable precision pipettes, preferably with disposable plastic tips. (A manifold multi-channel pipette is desirable for large assays.)
3. Cell Extraction Buffer (see Recommended Formulation, p. 10).
4. Distilled or deionized water.
5. Plate washer: automated or manual (squirt bottle, manifold dispenser, etc.).
6. Data analysis and graphing software. Graph paper: linear (Cartesian), log-log, or semi-log, as desired.
7. Glass or plastic tubes for diluting and aliquoting standard.
8. Absorbent paper towels.
9. Calibrated beakers and graduated cylinders in various sizes.

### **PROCEDURAL NOTES/LAB QUALITY CONTROL**

1. When not in use, kit components should be refrigerated. All reagents should be warmed to room temperature before use.
2. **Microtiter plates should be allowed to come to room temperature before opening the foil bags.** Once the desired number of strips has been removed, immediately reseal the bag and store at 2 to 8°C to maintain plate integrity.

3. Samples should be frozen if not analyzed shortly after collection. Avoid multiple freeze-thaw cycles of frozen samples. Thaw completely and mix well prior to analysis.
4. If particulate matter is present, centrifuge or filter prior to analysis.
5. All standards, controls and samples should be run in duplicate.
6. Samples that are greater than the highest standard point should be diluted with *Standard Diluent Buffer* and retested.
7. When pipetting reagents, maintain a consistent order of addition from well-to-well. This ensures equal incubation times for all wells.
8. Cover or cap all reagents when not in use.
9. **Do not mix or interchange different reagent lots from various kit lots.**
10. Do not use reagents after the kit expiration date.
11. Read absorbances within 2 hours of assay completion.
12. In-house controls should be run with every assay. If control values fall outside pre-established ranges, the accuracy of the assay is suspect.
13. All residual wash liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. ***Never*** insert absorbent paper directly into the wells.
14. Because *Stabilized Chromogen* is light sensitive, avoid prolonged exposure to light. Also avoid contact between *Stabilized Chromogen* and metal, or color may develop.



## SAFETY

All blood components and biological materials should be handled as potentially hazardous. Follow universal precautions as established by the Centers for Disease Control and Prevention and by the Occupational Safety and Health Administration when handling and disposing of infectious agents.

## DIRECTIONS FOR WASHING

**Incomplete washing will adversely affect the test outcome.** All washing must be performed with *Wash Buffer Concentrate (25X)* provided.

Washing can be performed manually as follows: completely aspirate the liquid from all wells by gently lowering an aspiration tip (aspiration device) into the bottom of each well. Take care not to scratch the inside of the well.

After aspiration, fill the wells with at least 0.4 mL of diluted wash solution. Let soak for 15 to 30 seconds, then aspirate the liquid. Repeat as directed under **ASSAY METHOD**. After the washing procedure, the plate is inverted and tapped dry on absorbent tissue.

Alternatively, the wash solution may be put into a squirt bottle. If a squirt bottle is used, flood the plate with wash buffer, completely filling all wells. After the washing procedure, the plate is inverted and tapped dry on absorbent tissue.

If using an automated washer, the operating instructions for washing equipment should be carefully followed. If your automated washer allows, 30 second soak cycles should be programmed into the wash cycle.

## PROCEDURE FOR EXTRACTION OF PROTEINS FROM CELLS

### A. Recommended Formulation of Cell Extraction Buffer:

10 mM Tris, pH 7.4

100 mM NaCl

1 mM EDTA

1 mM EGTA

1 mM NaF

20 mM  $\text{Na}_4\text{P}_2\text{O}_7$

2 mM  $\text{Na}_3\text{VO}_4$

1% Triton X-100

10% glycerol

0.1% SDS

0.5% deoxycholate

1 mM PMSF (stock is 0.3 M in DMSO)

Protease inhibitor cocktail (e.g., Sigma Cat. # P-2714) (reconstituted according to manufacturer's guideline). Add 100  $\mu\text{L}$  per 1 mL Cell Extraction Buffer.

This buffer is stable for 2 to 3 weeks at 4°C or for up to 6 months when aliquoted (without protease inhibitors and PMSF added) and stored at -20°C. When stored frozen, the Cell Extraction Buffer should be thawed on ice. This buffer (minus protease inhibitor cocktail and PMSF) can be obtained from Invitrogen Cat. # FNN0011. **Important:** add the protease inhibitors just before using. The stability of protease inhibitor supplemented Cell Extraction Buffer is 24 hours at 4°C. PMSF is very unstable and must be added prior to use, even if added previously.

## B. Protocol for Cell Extraction

This protocol has been successfully applied to several cell lines. Researchers should optimize the cell extraction procedures for their own applications.

1. Collect cells in PBS by centrifugation (non-adherent) or scraping from culture flasks (adherent).
2. Wash cells twice with cold PBS.
3. Remove and discard the supernatant and collect the cell pellet. (At this point the cell pellet can be frozen at -80°C and lysed at a later date).
4. Lyse the cell pellet in Cell Extraction Buffer for 30 minutes, on ice, with vortexing at 10 minute intervals. The volume of Cell Extraction Buffer depends on the cell number in cell pellet and expression of STAT5a [pY694]. For example,  $5 \times 10^7$  Jurkat cells grown in RPMI-1640 plus 10% FBS and treated with 100 ng/mL IFN- $\alpha$  for 15 minutes can be extracted in 1 mL of Extraction Buffer. Under these conditions, use of 1-10  $\mu$ L of the clarified cell extract diluted to a volume of 100  $\mu$ L/well in *Standard Diluent Buffer* (See **Assay Method**) is sufficient for the detection of STAT5a [pY694].
5. Transfer extract to microcentrifuge tubes and centrifuge at 13,000 rpm for 10 minutes at 4°C.
6. Aliquot the clear lysate to clean microfuge tubes. These samples are ready for assay. Lysates can be stored at -80°C. Avoid multiple freeze/thaw cycles.

## REAGENT PREPARATION AND STORAGE

### A. Reconstitution and Dilution of Hu STAT5a [pY694] Standard

**Note:** This *Hu STAT5a [pY694] Standard* is prepared from phosphorylated recombinant STAT5a C-terminal region. One Unit of standard is equivalent to 0.95  $\mu$ g of HEL cell lysate treated with 1 mM sodium vanadate for 6 hours.

1. Reconstitute *Hu STAT5a [pY694] Standard* with *Standard Diluent Buffer*. Refer to standard vial label for instructions. Swirl or mix gently and allow to sit for 10 minutes to ensure complete reconstitution. Label as 100 Units/mL STAT5a [pY694]. Use the standard within 1 hour of reconstitution.
2. Add 0.25 mL of *Standard Diluent Buffer* to each of 6 tubes labeled 50, 25, 12.5, 6.25, 3.12 and 1.6 Units/mL STAT5a [pY694].
3. Make serial dilutions of the standard as described in the following dilution table. Mix thoroughly between steps.

**B. Dilution of Hu STAT5a [pY694] Standard**

<b>Standard:</b>	<b>Add:</b>	<b>Into:</b>
100 Units/mL	Prepare as described in step 1	
50 Units/mL	0.25 mL of the 100 Units/mL std.	0.25 mL of the Diluent Buffer
25 Units/mL	0.25 mL of the 50 Units/mL std.	0.25 mL of the Diluent Buffer
12.5 Units/mL	0.25 mL of the 25 Units/mL std.	0.25 mL of the Diluent Buffer
6.25 Units/mL	0.25 mL of the 12.5 Units/mL std.	0.25 mL of the Diluent Buffer
3.12 Units/mL	0.25 mL of the 6.25 Units/mL std.	0.25 mL of the Diluent Buffer
1.6 Units/mL	0.25 mL of the 3.12 Units/mL std.	0.25 mL of the Diluent Buffer
0 Units/mL	0.25 mL of the Diluent Buffer	An empty tube

Remaining reconstituted standard should be discarded or frozen in aliquots at -80°C for further use. Standard can be frozen and thawed one time only without loss of immunoreactivity.

### C. Storage and Final Dilution of Anti-Rabbit IgG HRP (100X)

**Please Note:** The *Anti-Rabbit IgG HRP (100X)* is in 50% glycerol. This solution is viscous. To ensure accurate dilution, allow *Anti-Rabbit IgG HRP (100X)* to reach room temperature. Gently mix. Pipette *Anti-Rabbit IgG HRP (100X)* slowly. Remove excess concentrate solution from pipette tip by gently wiping with clean absorbent paper.

1. Dilute 10  $\mu\text{L}$  of this 100X concentrated solution with 1 mL of *HRP Diluent* for each 8-well strip used in the assay. Label as Anti-Rabbit IgG HRP Working Solution.

For Example:

# of 8-Well Strips	Volume of Anti-Rabbit IgG HRP (100X)	Volume of Diluent
2	20 $\mu\text{L}$ solution	2 mL
4	40 $\mu\text{L}$ solution	4 mL
6	60 $\mu\text{L}$ solution	6 mL
8	80 $\mu\text{L}$ solution	8 mL
10	100 $\mu\text{L}$ solution	10 mL
12	120 $\mu\text{L}$ solution	12 mL

2. Return the unused *Anti-Rabbit IgG HRP (100X)* to the refrigerator.

#### D. Dilution of Wash Buffer

Allow the *Wash Buffer Concentrate (25X)* to reach room temperature and mix to ensure that any precipitated salts have redissolved. Dilute 1 volume of the *Wash Buffer Concentrate (25X)* with 24 volumes of deionized water (e.g., 50 mL may be diluted up to 1.25 liters, 100 mL may be diluted up to 2.5 liters). Label as Working Wash Buffer.

Store both the concentrate and the Working Wash Buffer in the refrigerator. The diluted buffer should be used within 14 days.

#### ASSAY METHOD: PROCEDURE AND CALCULATIONS

**Be sure to read the *Procedural Notes/Lab Quality Control* section before carrying out the assay.**

Allow all reagents to reach room temperature before use. Gently mix all liquid reagents prior to use.

**Note:** A standard curve must be run with each assay.

1. Determine the number of 8-well strips needed for the assay. Insert these in the frame(s) for current use. (Re-bag extra strips and frame. Store these in the refrigerator for future use.)
2. Add 100  $\mu$ L of the *Standard Diluent Buffer* to zero wells. Well(s) reserved for chromogen blank should be left empty.
3. Add 100  $\mu$ L of standards, controls, and diluted samples (>1:10 dilution for cell extract) to the appropriate microtiter wells. Tap gently on side of plate to thoroughly mix. (See **REAGENT PREPARATION AND STORAGE**, Section B.)
4. Cover wells with *plate cover* and incubate for **2 hours at room temperature**.

5. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
6. Pipette 100  $\mu$ L of *Hu STAT5a [pY694] Detection Antibody* solution into each well except the chromogen blank(s). Tap gently on the side of the plate to mix.
7. Cover wells with *plate cover* and incubate for **1 hour at room temperature**.
8. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
9. Add 100  $\mu$ L *Anti-Rabbit IgG HRP Working Solution* to each well except the chromogen blank(s). (Prepare the working dilution as described in **REAGENT PREPARATION AND STORAGE**, Section C.)
10. Cover wells with the *plate cover* and incubate for **30 minutes at room temperature**.
11. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
12. Add 100  $\mu$ L of *Stabilized Chromogen* to each well. The liquid in the wells will begin to turn blue.
13. Incubate for **30 minutes at room temperature and in the dark**. **Please Note: Do not cover the plate with aluminum foil or metalized mylar.** The incubation time for chromogen substrate is often determined by the microtiter plate reader used. Many plate readers have the capacity to record a maximum optical density (O.D.) of 2.0. The O.D. values should be monitored and the substrate reaction stopped before the O.D. of the positive wells



exceed the limits of the instrument. The O.D. values at 450 nm can only be read after the *Stop Solution* has been added to each well. If using a reader that records only to 2.0 O.D., stopping the assay after 20 to 25 minutes is suggested.

14. Add 100  $\mu$ L of *Stop Solution* to each well. Tap side of plate gently to mix. The solution in the wells should change from blue to yellow.
15. Read the absorbance of each well at 450 nm having blanked the plate reader against a chromogen blank composed of 100  $\mu$ L each of *Stabilized Chromogen* and *Stop Solution*. Read the plate within 2 hours after adding the *Stop Solution*.
16. Plot on graph paper the absorbance of the standards against the standard concentration. (Optimally, the background absorbance may be subtracted from *all* data points, including standards, unknowns and controls, prior to plotting.) Draw the best smooth curve through these points to construct the standard curve. If using curve fitting software, the four parameter algorithm provides the best curve fit.
17. Read the STAT5a [pY694] concentrations for unknown samples and controls from the standard curve plotted in step 16. **Multiply value(s) obtained for sample(s) by the appropriate dilution factor to correct for the dilution with *Standard Diluent Buffer*.** (Samples producing signals higher than the highest standard [100 Units/mL] should be further diluted in *Standard Diluent Buffer* and reanalyzed, multiplying the concentration by the appropriate dilution factor.)

## TYPICAL DATA

The following data were obtained for the various standards over the range of 0 to 100 Units/mL STAT5a [pY694].

Standard STAT5a [pY694] (Units/mL)	Optical Density (450 nm)
100	3.15
50	1.77
25	1.18
12.5	0.78
6.25	0.44
3.12	0.29
1.6	0.24
0	0.15

## LIMITATIONS OF THE PROCEDURE

Do not extrapolate the standard curve beyond the 100 Units/mL standard point; the dose-response is non-linear in this region and accuracy is difficult to obtain. Dilute samples >100 Units/mL with *Standard Diluent Buffer*; reanalyze these and multiply results by the appropriate dilution factor.

The influence of various extraction buffers has not been thoroughly investigated. The rate of degradation of native STAT5a or dephosphorylation of STAT5a [pY694] in various matrices has not been investigated. Although STAT5 degradation or dephosphorylation of STAT5a [pY694] in the Cell Extraction Buffer described in this protocol has not been seen to date, the possibility of this occurrence cannot be excluded.

**For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.**


## PERFORMANCE CHARACTERISTICS

### SENSITIVITY

The analytical sensitivity of this assay is <1 Unit/mL of STAT5a [pY694]. This was determined by adding two standard deviations to the mean O.D. obtained when the zero standard was assayed 30 times.

The sensitivity of this ELISA was compared to Western blotting using known quantities of STAT5a [pY694]. The data presented in Figure 1 show that the sensitivity of the ELISA is approximately 2x greater than that of Western blotting. The bands shown in the Western blotting data were developed using rabbit anti-STAT5a [pY694], and an alkaline phosphatase conjugated anti-rabbit IgG followed by chemiluminescent substrate and autoradiography.

**Figure 1: Detection of STAT5a [pY694] by ELISA vs Western Blot: HEL Cells Stimulated by Sodium Vanadate**

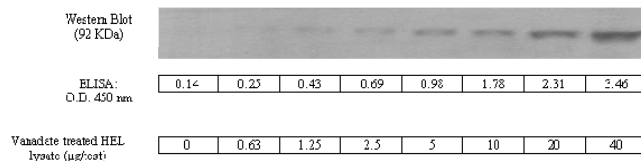

## PRECISION

### 1. Intra-Assay Precision

Samples of known STAT5a [pY694] concentrations were assayed in replicates of 12 to determine precision within an assay.

	Sample 1	Sample 2	Sample 3
Mean (Units/mL)	27.64	13.72	6.45
SD	1.66	1.35	0.57
%CV	6.01	9.84	8.83

SD = Standard Deviation

CV = Coefficient of Variation

### 2. Inter-Assay Precision

Samples were assayed 36 times in multiple assays to determine precision between assays.

	Sample 1	Sample 2	Sample 3
Mean (Units/mL)	27.49	13.72	6.45
SD	1.87	1.35	0.70
%CV	6.81	9.87	10.90

SD = Standard Deviation

CV = Coefficient of Variation


## RECOVERY

To evaluate recovery, STAT5a [pY694] Standard was spiked at 7 different concentrations into 10% Cell Extraction Buffer. The average recovery was 91%.

## PARALLELISM

Natural STAT5a [pY694] from sodium vanadate treated HEL cell lysate was serially diluted in *Standard Diluent Buffer*. The optical density of each dilution was plotted against the STAT5a [pY694] standard curve. Parallelism demonstrated by the figure below indicates that the standard accurately reflects STAT5a [pY694] content in samples.

**Figure 2**

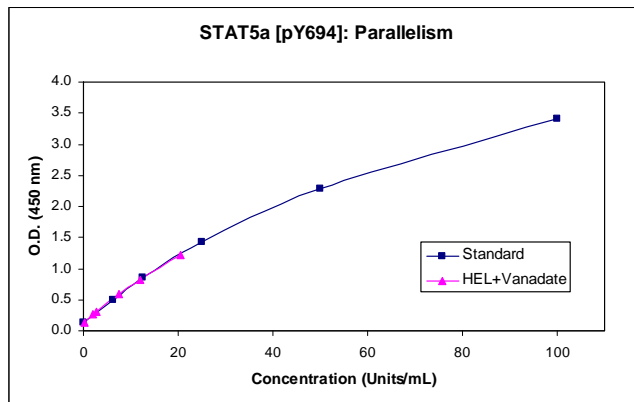

## LINEARITY OF DILUTION


HEL cells were grown in tissue culture medium containing 10% fetal bovine serum, treated with 1 mM sodium vanadate for 6 hours and lysed with Cell Extraction Buffer. This lysate was diluted in *Standard Diluent Buffer* over the range of the assay and measured for STAT5a [pY694]. Linear regression analysis of sample values versus the expected concentrations yielded a correlation coefficient of 0.99.

<b>Dilution</b>	<b>Cell Lysate</b>		
	<b>Measured (Units/mL)</b>	<b>Expected (Units/mL)</b>	<b>% Expected</b>
Neat	34.29	34.29	100
1/2	15.22	17.15	89
1/4	9.53	8.57	111
1/8	4.86	4.29	113
1/16	1.93	2.14	90

### SPECIFICITY

In Figure 3, HEL cell lysates were incubated with the capture antibody used in the STAT5a [pY694] ELISA (Cat. # KHO00761) (lane 2). An antibody specific for STAT5a and STAT5b was used as controls (lane 1 and 4, respectively). IgG beads were used as a negative control (lane 3). From these results the capture antibody only recognizes the a isoform of STAT5, but not the b isoform. Thus, the STAT5a [pY694] ELISA (Cat. # KHO00761) does not cross react with the STAT5b protein.


**Figure 3**

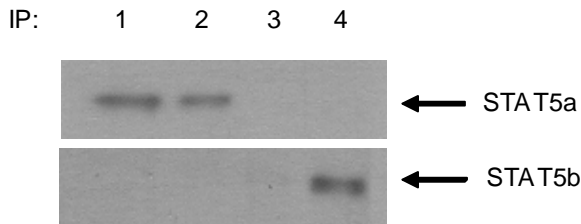




The specificity of this assay for phosphorylated STAT5a [pY694] was confirmed by peptide competition. The data presented in Figure 4 show that only the phospho-peptide containing the phosphorylated tyrosine 694 blocks the ELISA signal. The non-phosphorylated peptide sequence did not block the signal. Additionally, phosphorylated STAT1 peptide p701 did not block the signal, therefore confirming the specificity of the antibody for p694 only.

**Figure 4**

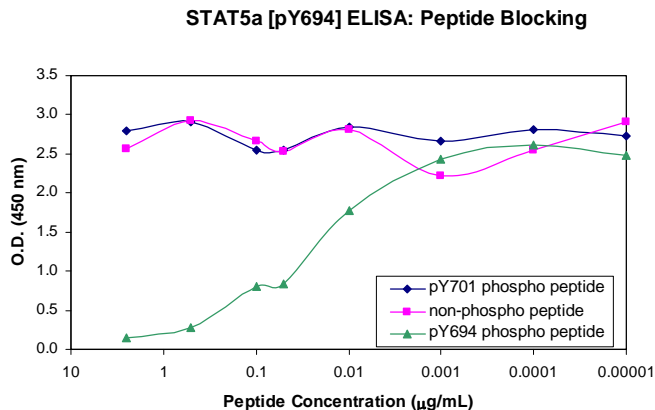

In Figure 5, HEL cells were treated with 1 mM sodium vanadate for 6 hours. Untreated cells were used as a negative control. Cell extracts were prepared and analyzed with the STAT5a [pY694] ELISA and Invitrogen's STAT5a (Total) ELISA (Cat. # KHO0751). The results show that the phosphorylation of STAT5a is increased in sodium vanadate treated HEL cells, whereas the total level of STAT5a remains relatively constant in treated vs. untreated control.

**Figure 5**

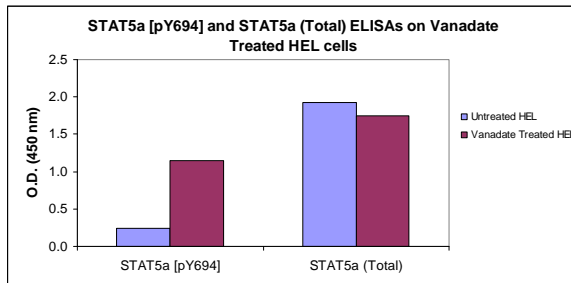














## REFERENCES

1. Tallquist, M., et al. (2004) PDGF signaling in cells and mice. *Cytokine Growth Factor Rev.* 15(4):205-13.
2. Gutierrez-Castellanos, S., et al. (2004) Differences in BCL-X(L) expression and STAT5 phosphorylation in chronic myeloid leukaemia patients. *Eur. J. Haematol.* 72(4):231-238.
3. Xu, W., et al. (2004) Geldanamycin, a heat shock protein 90-binding agent, disrupts Stat5 activation in IL-2-stimulated cells. *J. Cell. Physiol.* 198(2):188-196.
4. Rico-Bautista, E., et al. (2004) Downregulation of the growth hormone-induced Janus kinase 2/signal transducer and activator of transcription 5 signaling pathway requires an intact actin cytoskeleton. *Exp. Cell. Res.* 294(1):269-280.
5. Heath, C., et al. (2004) Critical role of STAT5 activation in transformation mediated by ZNF198-FGFR1. *J. Biol. Chem.* 279(8):6666-6673.
6. Schmitt-Ney M., et al. (1991)  $\beta$ -casein gene promoter activity is regulated by the hormone-mediated relief of transcriptional repression and a mammary-gland-specific nuclear factor. *Mol. Cell. Biol.* 11(7):3745-3755.
7. Wakao H., et al. (1992) Mammary gland-specific nuclear factor is present in lactating rodent and bovine mammary tissue and composed of a single polypeptide of 89 kDa. *J. Biol. Chem.* 267(23):16365-16370.
8. Wakao H., et al. (1994) Mammary gland factor (MGF) is a novel member of the cytokine regulated transcription factor gene family and confers the prolactin response. *EMBO J.* 13(18):2182-2191.

9. Welte T., et al. (1994) Prolactin-dependent activation of a tyrosine phosphorylated DNA binding factor in mouse mammary epithelial cells. *Mol. Endocrinol.* 8(8):1091-1102.
10. Shipley, J.M., et al. (2003) Down-regulation of STAT5b transcriptional activity by ligand-activated peroxisome proliferator-activated receptor (PPAR) alpha and PPARgamma. *Mol. Pharmacol.* 64(2):355-364.
11. Okutani, Y., et al. (2001) Src directly tyrosine-phosphorylates STAT5 on its activation site and is involved in erythropoietin-induced signaling pathway. *Oncogene* 20(45):6643-6650.
12. Gesbert, F., et al. (2000) Bcr/Abl activates transcription of the Bcl-X gene through STAT5. *Blood* 96(6):2269-2276.

**Important Licensing Information** - These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or our website, [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses. Unless otherwise indicated, these products are for research use only and are not intended for human or animal diagnostic, therapeutic or commercial use.

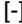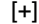**Explanation of symbols**

Symbol	Description	Symbol	Description
	Catalogue Number		Batch code
	Research Use Only		<i>In vitro</i> diagnostic medical device
	Use by		Temperature limitation
	Manufacturer		European Community authorised representative
	Without, does not contain		With, contains
	Protect from light		Consult accompanying documents
	Directs the user to consult instructions for use (IFU), accompanying the product.		

Copyright © Invitrogen Corporation. 24 February 2010





## STAT5a [pY694] Assay Summary

Incubate 100  $\mu$ L Standard or Cell Extract (>1:10)  
for 2 hours at RT



**aspirate and wash 4x**

Incubate 100  $\mu$ L of Detection Antibody  
for 1 hour at RT



**aspirate and wash 4x**

Incubate 100  $\mu$ L of HRP Anti-Rabbit Antibody  
for 30 minutes at RT



**aspirate and wash 4x**

Incubate 100  $\mu$ L of Stabilized Chromogen  
for 30 minutes at RT



Add 100  $\mu$ L of Stop Solution and read at 450 nm



**Total time: 4 hours**

