


Superior cloning technologies

Directional TOPO® Cloning

The most effective technology available for cloning DNA

In the 5-minute Directional TOPO® Cloning reaction, the enzyme topoisomerase I is used in place of DNA ligase, greatly increasing the success and speed of the cloning reaction. Directional TOPO® Cloning allows the cloning of PCR products in a specific 5' to 3' orientation.

Directional TOPO® Cloning procedure

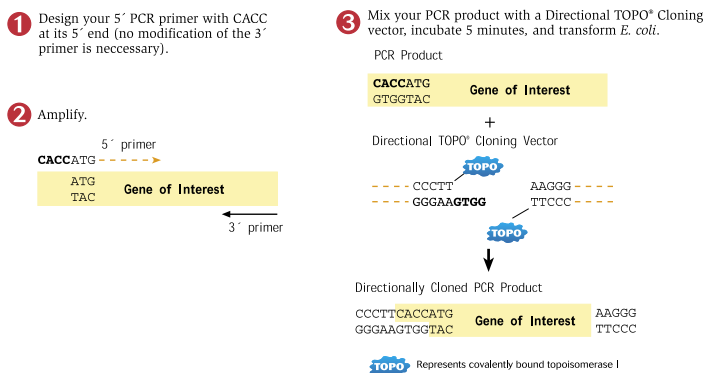


Gateway® Cloning Technology


The universal cloning and expression platform

Get rapid, highly efficient cloning of DNA segments across multiple systems and vectors with Gateway® Technology. After generating an entry clone, you can quickly and easily move a gene into different expression vectors.

The flexibility of Gateway® Technology

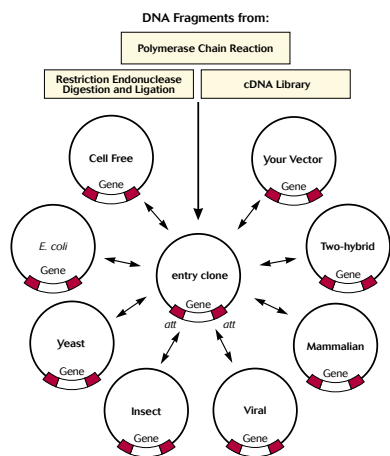

The Gateway® LR recombination reaction

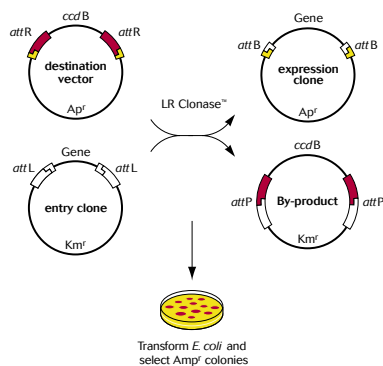


Advanced labeling and detection tool

Lumio™ Technology

Specific, sensitive, and rapid detection of protein expression

The versatile Lumio™ platform uses a small, six-amino acid novel recognition sequence and target-specific fluorescent labeling reagents, enabling you to quickly detect proteins in gels using a standard UV transilluminator or laser based scanner. It eliminates protein staining and time-consuming western blot procedures.

Lumio™ detection of CAT under UV light (left) and Coomassie® stained (right)

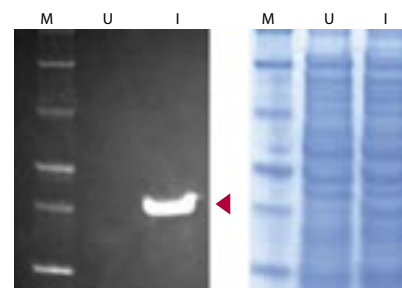

M: Fluorescent marker
U: pET161-DEST/CAT uninduced
I: pET161-DEST/CAT Induced

CAT was expressed from the Champion™ pET161-DEST vector. Samples were run on a 12% NuPAGE® Bis-Tris gel with MES running buffer with the Lumio™ Green Detection Reagent and optimized Lumio™ Gel Sample Buffer. The gel was visualized under ultraviolet light (left) and then stained with Coomassie® Blue (right).

Diverse Champion™ pETvector options for *E. coli* expression

Vector	Cat. no.	Position	Tag	Cleavage Protease	Antibiotic resistance	Benefit
pET100/D-TOPO®	K100-01	N-term	6xHis-Xpress™	EK	Amp	Cleavable detection and purification tag
pET101/D-TOPO®†	K101-01	C-term	V5-6xHis	-	Amp	Detection and purification tag
pET102/D-TOPO®	K102-01	N-term	Thioredoxin	EK	Amp	Cleavable thioredoxin tag enhances protein translation and solubility
		C-term	V5-6xHis	-	-	Detection and purification tag
pET151/D-TOPO®	K151-01	N-term	6xHis-V5	TEV	Amp	Cleavable thioredoxin tag enhances protein translation and solubility
pET160/GW/D-TOPO®	K160-01	N-term	6xHis-Lumio™	TEV	Amp	Direct fluorescent in-gel detection and cleavable purification tag
pET161/GW/D-TOPO®	K161-01	C-term	Lumio™-6xHis	-	Amp	Direct fluorescent in-gel detection and cleavable purification tag
pET200/D-TOPO®	K200-01	N-term	6xHis-Xpress™	EK	Kan	Cleavable detection and purification tag
pET SUMO	K300-01	N-term	6xHis-SUMO	-	Kan	Highest protein solubility with the SUMO tag; cleavable detection and purification tag; generation of native proteins and peptides
pET-DEST42	12276-010	C-term	V5-6xHis	-	Amp	Detection and purification tag
pET160-DEST	12583-035	N-term	6xHis-Lumio™	TEV	Amp	Direct fluorescent in-gel detection and cleavable purification tag
pET161-DEST	12583-043	C-term	Lumio™-6xHis	-	Amp	Direct fluorescent in-gel detection and purification tag
pET104-DEST	K104-01	N-term	BioEase™	EK	Amp	<i>in vivo</i> biotinylation tag for detection and purification with streptavidin

† Useful for native protein expression by inserting a stop codon before the C-terminal fusion sequence.

Improved expression host strains

Recombinant protein expression strains

Product	Cat. no.	Benefits
BL21 Star™(DE3)	C6010-03	Maximize protein production in BL21 <i>E. coli</i> by enhancing mRNA stability (see β-gal expression compared in BL21 and BL21 Star™ <i>E. coli</i> systems).
BL21 Star™(DE3)pLysS	C6020-03	Maximize protein production in BL21 <i>E. coli</i> . Addition of pLysS lowers basal-level expression for expression of proteins that are slightly growth inhibitive to <i>E. coli</i> .
BL21-AI™	C6070-03	Tight regulation and strong expression of proteins from the T7 promoter. Recommended for expression of toxic genes.

Efficient removal of fusion tags

AcTEV™ Protease: a stabilized TEV protease

Recognition Sequence		
Glu Asn Leu Tyr Phe Gln Gly ^TEV/AcTEV cleavage site		
Features	Quantity	Cat. no.
• Highly specific cleavage	1,000 units	12575-015
• Active from 4°C to 30°C and from pH 6.0 to pH 8.5	10,000 units	12575-023
• Increased enzyme stability for prolonged activity		
• 6xHis sequence facilitates removal following cleavage		

SUMO (ubiquitin-like protein) Protease

Features	Quantity	Cat. no.
• Highly specific cleavage activity*	100 units	12588-018
• Recognizes tertiary structure of SUMO		
• Active from 2°C to 37°C		
• 6xHis sequence facilitates removal following cleavage		
• Does not leave extra amino acids from protease recognition sequence following cleavage		

*The SUMO Protease can cleave any fusion to the SUMO protein fusion, except fusions that begin with proline.

EKMax™ Enterokinase, recombinant

Recognition Sequence		
Asp Asp Asp Asp Lys ^EK cleavage site		
Features	Quantity	Cat. no.
• Highly specific cleavage activity	250 units	E180-01
• Does not leave extra amino acids from protease recognition sequence following cleavage	1,000 units	E180-02
• Uses EK-Away™ Resin for protease removal		

Convenient media and selection

imMedia™ pre-mixed, pre-sterilized media: complete media in 5 minutes (no autoclaving of water or glassware required)

Product	Quantity	Cat. no.
For the preparation of liquid medium*		
imMedia™ Ampicillin Liquid	20 pouches	Q600-20
imMedia™ Kanamycin Liquid	20 pouches	Q610-20
imMedia™ Zeocin™ Liquid	20 pouches	Q620-20
For the preparation of agar plates**		
imMedia™ Ampicillin Agar	20 pouches	Q601-20
imMedia™ Kanamycin Agar	20 pouches	Q611-20
imMedia™ Zeocin™ Agar	20 pouches	Q621-20
For the preparation of agar plates with IPTG and X-gal**		
imMedia™ Ampicillin Blue	20 pouches	Q602-20
imMedia™ Kanamycin Blue	20 pouches	Q612-20

Effective concentrations of imMedia™ components: ampicillin (100 µg/ml), kanamycin (50 µg/ml), Zeocin™ (25 µg/ml), IPTG (100 µg/ml), and X-gal (100 µg/ml).

*Each pouch contains sufficient reagents to prepare 200 ml of media.

**Each pouch contains sufficient reagents to prepare 8-10 100 mm agar plates.

Traditional media

Product	Quantity	Cat. no.
LB Agar (Lennox L Agar)	500 g	22700-025
LB Broth (1X) (liquid)	500 ml	10855-021
LB Broth Base (Lennox Broth Base)	500 g	12780-052
Luria Broth Base (Miller's LB Broth Base)	500 g	12795-027
SELECT Agar	500 g	30391-023
S.O.C. Medium (liquid)	10 X 10 ml	15544-034
Terrific Broth	500 g	22711-022

Media components

Product	Quantity	Cat. no.
X-gal	1 g	15520-018
Bluo-gal (produces darker blue than X-gal)	1 g	15519-028
IPTG	1 g	15529-019

Antibiotics

Product	Quantity	Cat. no.
Ampicillin	200 mg	11593-019
Carbenicillin, disodium salt	5 g	10177-012
Kanamycin sulfate	5 g	11815-024
Kanamycin sulfate (liquid) (100X)	100 ml	15160-054
Streptomycin sulfate	100 g	11860-038
Zeocin™	1 g	R250-01

For more additional products and detailed information, visit www.invitrogen.com or call a Technical Service representative at 800 955 6288, option 2.

Zeocin is a trademark of CAYLA. These products may be covered by one or more Limited Use Label Licenses (See Invitrogen Catalog or www.invitrogen.com). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

©2004 Invitrogen Corporation. All rights reserved. Reproduction forbidden without permission. Printed in the U.S.A.



www.invitrogen.com

B-13508-r1/US 1104 x0xM